EDITO

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We are pleased to introduce our new catalogue.

Thanks to the close relationship that we maintain with our customers and partners, we were able to develop a wide range of products for many applications. The pages which follow contain more than a thousand references and a lot of information, including details about the proper use of devices.

More than a catalog, this book provides guidelines throughout your analyses and measurements. Designed to make your research easier when choosing your laboratory equipment, this handbook brings together our entire range. You will find:

- the Dujardin-Salleron's equipment – stills, ebulliometers, titration devices... a mix of elegance and functionality;
- the timeless instruments that became essential: mustimeters, alcoholometers, wine thieves... Their reliability and long-lasting quality never stop seducing;
- the necessary articles for your day-to-day analyses;
- a selection of products from renowned manufacturers along our own production;
- services of technical training, calibration, maintenance...

The « Made in France » by Dujardin-Salleron remains the must-have, synonymous with high-tech and sustainability. Let us say it another way: you can rely upon our products.

Committed to a process of continuous improvement, we wish to offer you a high quality of support and assistance. We hope you find this support useful. Your comments and suggestions are welcome to further improve our quality of service.

The entire team of Dujardin-Salleron Laboratories

The precursors of oenological analyses were Baumé, Gay-Lussac, and then came Pasteur. In the 18th and 19th centuries, they prepared the theoretical bases of fermentation. Following their famous wake, Jules Salleron became the forerunner in precision instruments, inventor of the first still, the acetimeter and above all the ebulliometer, never equalled till today. Jules Dujardin joined him very quickly, and the tandem kept on innovating for the useful engineering field. It brought to the Dujardin-Salleron label a scientific aura, that their successors have always been very keen on perpetuating. The company is not a family business any more since 1987, and this is to even better excel in the precision instrumentation field, serving œnology and spirits.

Including a multilingual sales and technical team able to answer to all your requests, a R&D department at the cutting edge of innovation, all our teams work to offer you the best testing instrumentation. To satisfy the very specific needs of professionals, we always bring appropriate technical and commercial solutions, even if they need to be invented or created. The renewal of ISO 9001:2015 certification is a supplementary guarantee of our commitment to provide quality goods and services to our partners.

Now assisted by a metrology laboratory named E.R.IN, we strengthen our expertise in the calibration, control, verification and maintenance of your instruments. It is essential for us to ensure the control of the instruments and offer you a high quality service throughout their whole lifetime.

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Tel : +33(2) 47 25 58 25 - info@dujardin-salleron.com - www.dujardin-salleron.com

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SERVICES D'ETALONNAGE CALIBRATION SERVICES

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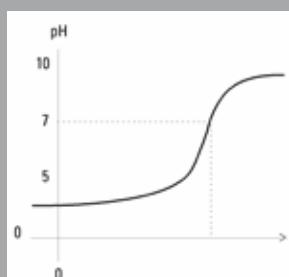
pH / Titration

pH is a crucial parameter in oenology, as it is a good indicator for acidity. It helps to adjust the harvest date and to know the corrections to make during vinification process.

Principle

A solution is systematically defined by its potential hydrogen (pH). It measures the chemical activity of protons (or H^+ hydrogen ions) in solution. It is often associated to acidity or basicity of a solution. In an aqueous solution at 25°C :

- A solution with $pH < 7$ is defined as acidic ; the lower its pH, the more acidic the solution.
- A solution with $pH = 7$ is defined as neutral (water is at $pH = 7$) ;
- A solution with $pH > 7$ is defined as basic ; the higher its pH, the more basic the solution.

Application

Titrimetry is a method used in analytical chemistry to determine the concentration of a chemical in solution. The titration is the implementation of this determination method, involving a chemical reaction between the chemical to measure and a standard reagent that is added progressively. Reaction must be total and measurable (pH, conductivity, color...).

pH metric titrations occur between acids and bases, with an end point detected by pH measurement or color indicator. All acids and bases can be titrated by standard solutions of strong bases or acids.

- Acidimetric titration : the user drops a base into an acidic volume. Total acidity in musts and wine, which mainly consists of tartaric acid and malic acid, is determined with this method. A standard solution of sodium hydroxide is dropped into a must or wine sample to bring pH to 7. The end point is determined by a color indicator (bromothymol blue) or by a pH meter (Titra EVO). The volume of sodium hydroxide used allows to calculate the acidity content of the sample.
- Alkalimetric titration : the user drops an acid into a basic sample. This method is not used in oenology.

Titration by redox potential involves another chemical reaction (more details p.14).

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Kit for determination of total acidity. It includes a volumetric graduated flask, a reagent and a standard solution, and a filtering set for musts. Fast and simple operations

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Complete kit for total acidity titration. It includes a 25 ml burette, a stand, a pipette, a crystallizing dish, a reagent and a standard solution for a direct readout.

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Complete kit in plastic case for total acidity titration. It includes an automatic 15 ml burette, a stand, a pipette, a crystallizing dish, a reagent and a standard solution in flexible flask for a direct readout.

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Dimensions (HxLxP), poids : 670 x 185 x 230 mm, 1,9 kg

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TA01 titration unit is designed to make easier the determinations by adding a titration solution with a changing colored indicator. The unit is composed of a 10 ml automatic burette in 1/20ml, controlled by an electrovalve, a magnetic stirrer with variable speed, and a retrolighted tray which offers a better visualization of the end of the titration.

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Descriptif complet p.16

TITRA EVO. This device enables pH measurement, as well as total acidity determination and volatile acidity after extraction. It avoids the difficult estimation of the detection of color change at the end point, and offers optimum accuracy and reproducibility of the results.

Supplied with accessories and products for around a hundred analysis.

Characteristics :

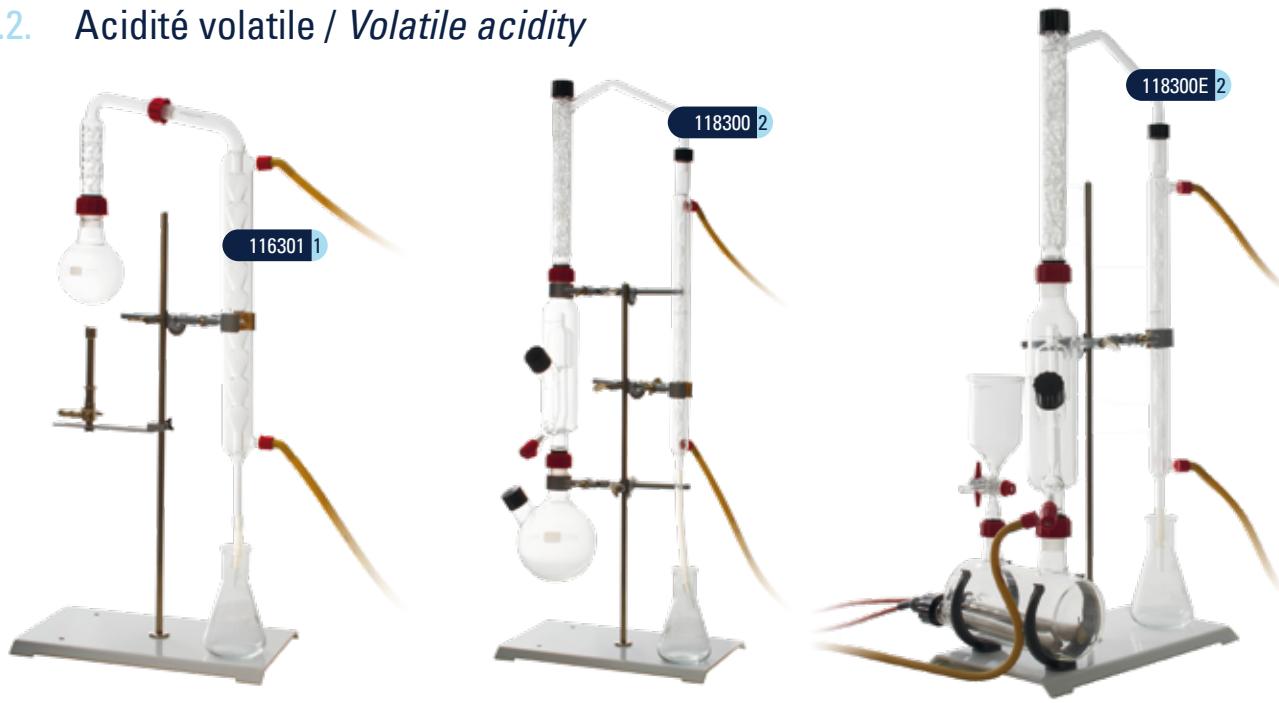
- Automatic titration
- Optional sample changer 15 positions
- Instinctive touch screen 7"
- Syringe 20 ml
- Memory 4 Go
- Data transfer on flash drive
- Possibility to combine to a feature of titration of SO₂

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Apparatus for the extraction of volatile acidity by direct distillation (Duclaux-Gayon method). Laboratory apparatus, with pyrex glassware, simple assembly/dismantling thanks to Rodavis system. Equipped with a rectifying column to stop lactic acid vapors.

116300 Appareil livré sans chauffage

Apparatus delivered without heating accessories

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The following apparatus are designed for the extraction of volatile acidity in conformity with European regulations (Official Journal L272 of 03/10/90) and the O.I.V standards such as those described in the Recueil des Méthodes Internationales d'Analyse des Vins et des Moûts.(Compendium of International Methods of Analysis of Wines and Musts).

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Apparatus for the extraction of volatile acidity by steam distillation (Mothe type). This apparatus with stand is composed of a boiler, an extraction flask, a rectifying column to stop lactic acid vapors and a cooling coil. Extraction time : around 8 minutes

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Automatic extractor Vola 2000. This apparatus is equipped with a system of instantaneous production of steam (without boiler under pressure) in highly resistant machine-cut aluminium. The high built-in electric heater (2800 W) enables the production of steam by injection, thanks to a dosage pump, of the water necessary for the generator throughout the entire extraction process.

Quick to start and reliable results make the VOLA 2000 advisable for laboratories dealing with wine analysis, spirits, fermented beverages, fruit juices and numerous other agriculture and food industry products (tomato sauces, vinegars, etc...).

The instrument is delivered with all necessary glassware and products for the extraction (except demineralized water).

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Complete kit for raw volatile acidity titration. It includes a 25 ml burette, a stand, a pipette, a crystallizer, a reagent and a standard solution for a direct readout.

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Complete kit in plastic case for raw volatile acidity titration. It includes an automatic 15 ml burette, a stand, a pipette, a crystallizer, a reagent and a standard solution in flexible flask for a direct readout.

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Complete kit for raw and corrected volatile acidity, including two 25 ml burettes, a stand, a pipette, a crystallizer, an Erlenmeyer flask, standard solutions and reagents.

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Complete kit for acetic acidity titration in vinegars. It includes a graduated burette in acetimetric degrees, a stand, a pipette, a crystallizer, a reagent and a standard solution for a direct readout.

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Complete kit in plastic case for acetic acidity titration in vinegars. It includes an automatic 15 ml burette, a stand, a pipette, a crystallizer, a reagent and a standard solution in flexible flask for a direct readout.

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Complete kit for chromatography on paper. It includes an amber glass ground joint flask, Whatman chromatographic paper, micro-pipette, reagents and standard samples.

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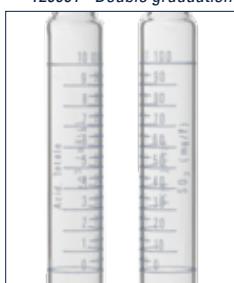
120601 - Double graduation
120601 - Double graduation

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SO₂ and total acidity kit. It enables to quickly estimate free SO₂ and total acidity. It contains a volumetric graduated flask, a reagent and a standard solution for each titration. Stable solutions, fast and simple operations.

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Complete kit for free and total SO₂ titration (Ripper method). It includes a 25 ml burette, a stand, pipettes, an Erlenmeyer flask, reagents and a standard solution.

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Equipped with an automatic burette, engraved and graduated in mg/l of SO₂, on amber glass flask, with rubber bulb and siphon, for a direct readout.

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Complete kit in plastic case for the titration of free and total SO₂. Free and total SO₂ are successively titrated by a stable standard solution of Iodide/Iodate (KIKIO₃) , which progressively gives iodine during titration. All volumes required can be measured with the specific measuring cylinder.

The case contains an automatic 25 ml burette, a stand, a measuring cylinder, an Erlenmeyer flask, reagents and a standard solution in flexible flask.

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SO₂ determination

SO₂, or sulfurous anhydride, is widely used in oenology. At any step of winemaking process, in particular at the end of alcoholic and malo-lactic fermentations, during ageing and before bottling, it has to be controlled. Its determination is done by titration, directly or after extraction.

By titration

This titration is based on a redox reaction, during which a transfer of electrons occurs. The chemical that loses an electron is called "reducteur" and the one that wins one is called "oxydant". Iodine is used as an oxidizing solution to determine SO₂ in acidic conditions.

After all dissolved SO₂ have reacted, iodine is in excess in the solution. This end point is detected by :

- Colorimetry, or Ripper method : a starch solution indicates the end of titration. Added to the sample before the titration, it changes from colorless to blue in the presence of iodine.

- Potentiometry : a double platinum electrode detects precisely the end of titration (Sulfilyser+ and TITRA EVO).

In both cases, the volume of iodine used is multiplied by a coefficient, automatically or by the user, to know the content of SO₂ (cf. chart p.121). Free SO₂ is determined directly, while total SO₂ is determined after releasing combinations of SO₂ with a concentrated solution of sodium hydroxide.

By aspiration-oxidation

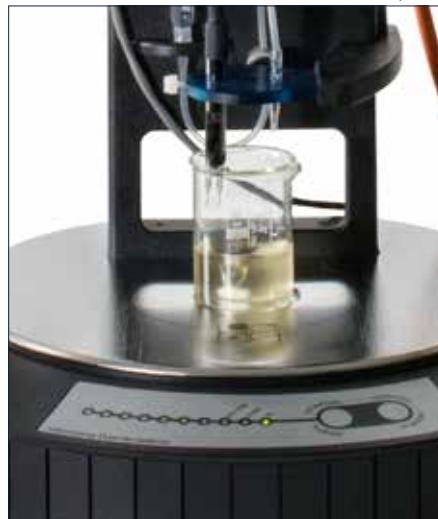
The reference method of OIV/EC recommends the use of a Frantz Paul apparatus, characterized by specific dimensions. SO₂ is distilled in acidic conditions by an extraction with gas flow (air or nitrogen). It is collected into an oxidizing solution of hydrogen peroxide (H₂O₂) in which it is converted into sulfuric acid (H₂SO₄), and then quantified by acidimetry.

Free SO₂ is extracted without heating, while total SO₂ is extracted after applying heat to the sample, to enable the extraction of combined SO₂. This method has the advantage of avoiding the influence of the various compounds of the sample, but it requires a rigorous laboratory practice.

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Sulfilyser+. Semi-automatic titrator to determine easily and quickly free and total SO2. It avoids the difficult estimation of the end of titration with Ripper method.

The Sulfilyser + is particularly adapted to highly colored red wines, lies for distillation, fruit juices, alcoholic drinks and vinegars. After the preparation of the sample, the titration is automatic until all the LED turn on, showing the end of the titration. The result is obtained multiplying by a coefficient the volume of solution used to be read on the burette.

The instruments is supplied with the titration glassware and the chemicals needed for approximately 100 analysis.

Characteristics :

- Automatic titration
- Excellent repeatability
- Integrated magnetic stirrer
- Removable stainless steel tray to make the cleaning easier
- Linear guidance of the support of the agitator and the probe
- Auto-zeroing 10 ml burette, graduated 1/20
- Power supply : 110-220V - 50/60 Hz – CE
- Dimensions (HxWxD), weight : 510 x 240 x 350 mm, 2,7 kg

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TITRA EVO. Automatic multiparameter titrator for quick determination of free and total SO₂, pH, volatile acidity on distillate. The device automatically adds the reagents before the titration. It is available with an autosampler to analyse up to 15 samples in a row. Its intuitive touch screen makes it particularly easy to use.

Iodometric determination of SO₂

With a double platinum electrode that detects an electric current in the sample, the instrument performs the titration and stops automatically when the oxidizing solution is in excess. The device calculates the result depending on the selected programme (free or total SO₂) and displays it directly after a few seconds.

Determination of pH and acidity

This device allows to determine pH, total acidity directly in the sample, and volatile acidity after extraction, without any need to see a color change at the end of titration. This enables optimum accuracy and reproducibility of the results.

Supplied with titration glassware and the chemicals needed for over a hundred analyses.

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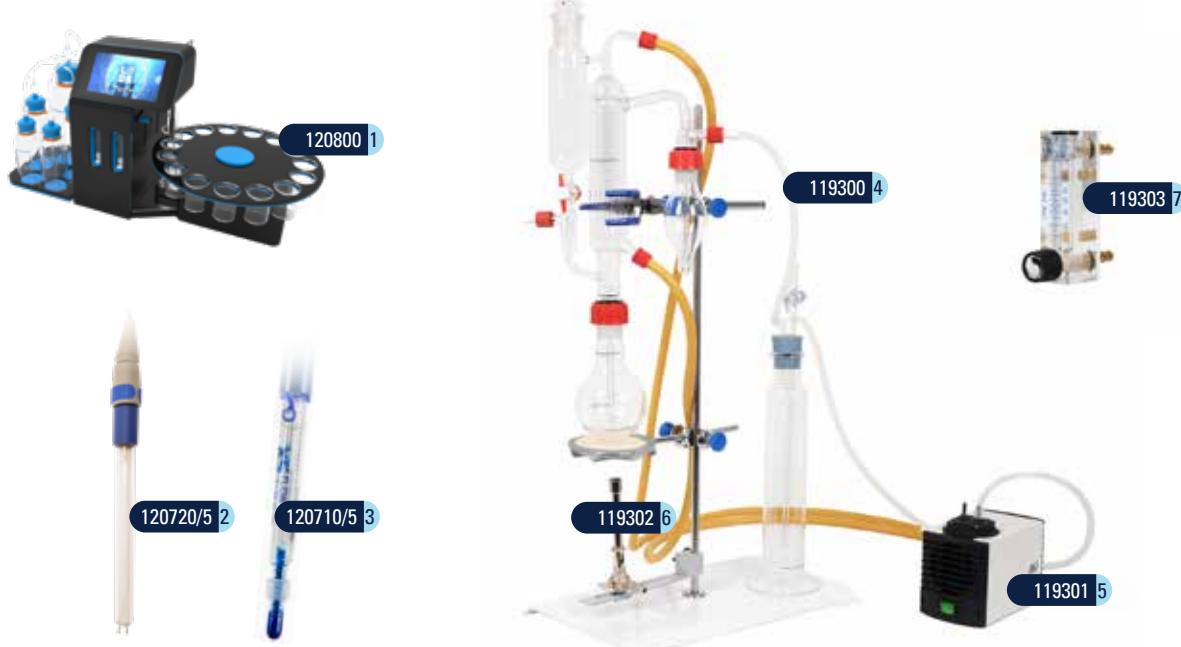
* possible evolution after initial purchase

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Sample changer 15 positions. Supplied separately, it is easy to connect to Titra EVO, to place and to activate, even on a device already set-up.

- Automatic rinsing of the electrode
- Integrated magnetic stirring
- Time per sample : free SO₂ approx.30s, total acidity approx.60s
- Dimensions Titra EVO + changer : 75 x 36 x 44 cm

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2 120720/5 Electrode double de platine (câble non inclus)
Double platinum electrode (without cable)

3 120710/5 Electrode pH (câble non inclus)
pH electrode (without cable)

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Frantz Paul apparatus. Apparatus for the extraction of free and total SO₂ by aspiration-oxydation according to the reference method OIV/CE. The apparatus is composed of a stand, a ground joint extraction flask, a cooler unit, a ground joint bubbling flask and a cylinder. Supplied without heating nor chemicals.

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Electric vacuum mini-compressor. Supplied with tube and adapters for Frantz Paul.

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Heating and fittings for Frantz Paul apparatus. Includes a gas burner, a ring, a gauze and holders.

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Flowmeter for Frantz Paul. 0 to 1 litre/minute. Delivered with connectors.

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FP auto. Automatic Frantz Paul apparatus for the fast extraction of free and total SO₂ by aspiration-oxydation according to the reference method OIV/CE.

Characteristics :

- Selection of the operating time with a timer
- Automatic stop combined to the closing of the circulating cooling system
- Built-in heating mantle for the extraction of total SO₂
- Vacuum pump integrated to the apparatus
- Adjustable air flow thanks to an built-in flowmeter
- Dimensions (HxWxD) : 70 x 37 x 34 cm

Supplied with glassware and chemicals needed for the SO₂ extraction. Available in 110V upon request.

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Titration kit for Frantz Paul. It enables to directly titrate the sulfuric acid collected in the bubbling flask, with a freshly diluted solution of sodium hydroxide. This kit contains a 10 ml burette, a stand, a standard solution and the accessories for the dilution.

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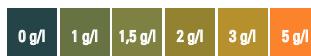
Fermentest®. Simple and quick method to determine reducing sugars in must and wine. It quickly and easily differentiates dry wines from wines with higher sugar content, by comparison of the color obtained after reaction to a color chart.

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Complete kit for determination of residual sugars in wines and musts with Fehling method. It contains a burette on a stand, a heating system, products for decoloration, Fehling liquor, Erlenmeyer flask and a wooden clamp.

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Ebulliometry

Studied and developed by the founders of Laboratoires Dujardin-Salleron, the ebulliometry is a simple and direct method to determine the alcohol content of dry wines and vinegars. It is used worldwide not only by winemakers, but also in other fields of food industry.

Principle

The boiling point is one of the characteristics of a liquid : at a pressure of 1 atmosphere, water boils at 100°C and ethanol at 78,4°C. As wine is a solution mainly composed of water and alcohol, its boiling point depends on its alcohol content. The lower its boiling point, the higher its alcohol content.

Atmospheric pressure and altitude influence the boiling point. Therefore, a calibration is required to determine the environment conditions at the moment of the measurement. So the first step is to measure the boiling point of water, or of a standard wine, to adjust the correspondence table "boiling point / alcohol content". Then the boiling point of the sample can be measured and reported to the correspondence table to know the alcoholic degree.

Sugars in solution have an influence on boiling point. Below 2 g/l, this influence is not important enough to reduce accuracy. Above this threshold, a correction must be applied, or another method must be used.

Recommendations

The measurement of boiling point requires the use of an ebulliometer, traditional or electric. In all cases, a measurement (excluding calibration) lasts about 7-8 minutes. This method is valid for dry wines below 20% A.b.V. with an accuracy of approximately $\pm 0,1\%$ A.b.V.

To obtain more accurate results, some precautions must be taken :

- use the same thermometer or the same probe to perform calibration and measurement ;
- respect scrupulously the volumes of standard solution and sample ;
- work with liquids and an ebulliometer at room temperature ;
- check that refrigeration is efficient to avoid loss of alcohol vapors during heating ;
- make sure there is no sugar in the sample ;
- renew calibration on a regular basis if the ebulliometer is not equipped with a sensor of atmospheric pressure.

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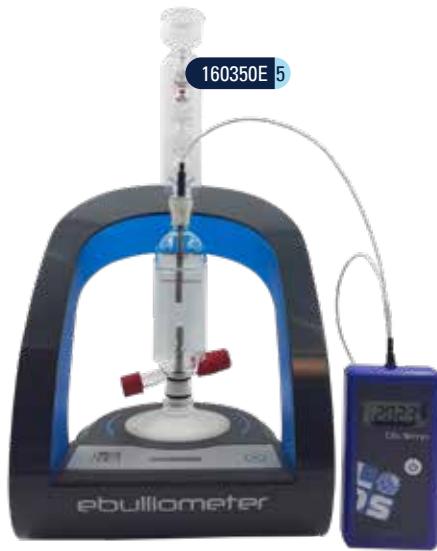
Complete ebulliometer in wooden case, alcohol heating.

Developed around 1870, continuously improved and recognised throughout the world, the Dujardin-Salleron Ebulliometer is the simplest instrument for directly and accurately (0,1% AbV) testing the alcoholic content of dry wines with normal composition.

Comes complete with wooden case, DS Temp precision thermometer, calculation disk, alcohol lamp, water/wine measurement tube and a box of wicks.

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Special ebulliometer for vinegar testing.

The same, with a boiler, the inner surface of which is covered with a protective coating tin/nickel.

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3 690114 Tube eau-vin/ Water-wine measurement tube

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Ebulliometer with DS Temp. With an electronically regulated electric heater and continuous refrigeration, the Ebulliometer enables to accurately determine the alcohol content of dry wines in about 7 minutes. With an ergonomic and original conception, the device is easy to set up and use. It is supplied with all needed accessories to make measurements immediately.

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Ebulliometer. With an electronically regulated electric heater and continuous refrigeration, the Ebulliometer enables to accurately determine the alcohol content of dry wines in about 7 minutes.

Thanks to the automatic atmospheric pressure compensation, there is no need of frequent calibrations.

With an ergonomic and original conception, the device is easy to set up and use. It is supplied with all needed accessories to make measurements immediately.

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Additional methods containing instruments and calculation formula to correct the result of the Ebulliometer according to the density of the sample.

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Ebullilog. Calculation software to obtain the Alcohol by Volume according to the boiling point and calibration.

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Ebulliotronic. With the same design as the Ebulliometer, this device is equipped with a boiler, the shape and dimensions of which are optimised. Thanks to the integrated calculation tables, the alcohol content is directly displayed in % AbV.

Sample types : dry wines, vinegars, kombucha.

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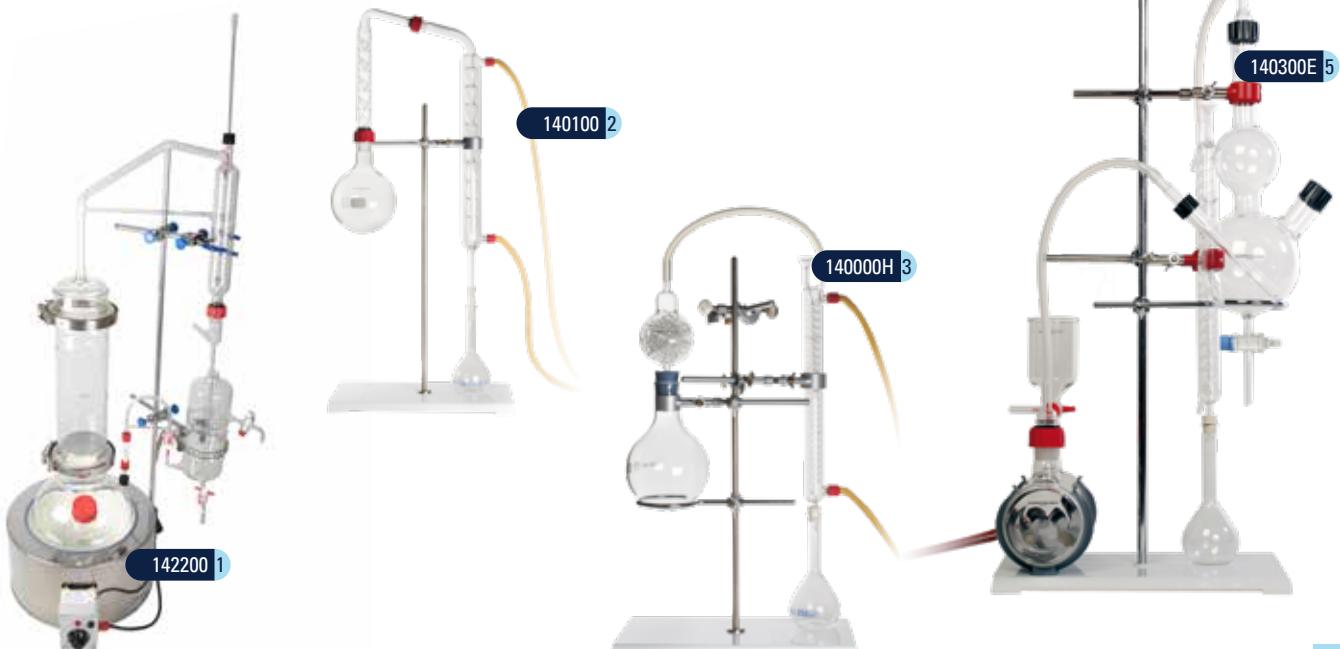
Still for plants and liquors. Glass apparatus to make tests by steam distillation of plants and infusions, conceived for distillers and perfumers. It works with a heating mantle, not provided (contact us).

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Dujardin-Salleron laboratory still. The Rodaviss system allows a quick and easy dismantling and makes the apparatus airtight. Equipped with a rectifying column and a reflux condenser.

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Mothe laboratory still for all degrees. Equipped with a rectifying bowl, a Graham condenser, and a stopper which ensure a perfect airtightness between the collecting flask and the condenser, it is specially adapted to the analysis of spirits over 25% AbV. Heating parts not included, to be ordered separately.

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Mothe distiller-extractor. It allows a quick steam extraction of alcohol and volatile acids, according to the official method. It is adapted to any kind of wine, liquor (<20% AbV), musts, even if there are sugars or solid pieces.

Equipped with a stand and a Graham condenser. Delivered without heating accessories. It is recommended to use a high flow gas burner Ø 40 mm.

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Mothe distiller-extractor with electric heating. Same principle and same applications as the apparatus 140300. Delivered with electric heating 230V - 1000W.

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Distillation - Extraction

Fractional distillation is the type I method recommended by OIV to obtain the Alcohol by Volume (A.b.V.) and volatile acidity, and type IV method for the sorbic acid. When determining the A.b.V., this method is more demanding than ebulliometry, but it is more accurate and it avoids the influence of other compounds on the sample, whatever the sugar and alcohol content.

Principle of laboratory distillation

Distillation is a process used to separate the various liquids in solution : they evaporate one after another, and their vapor is condensed in a condenser to collect a distillate. Laboratory distillation is called "fractional" because of the insertion of a rectifying column in order to achieve a better separation of the liquids of the mixture during extraction.

Applications

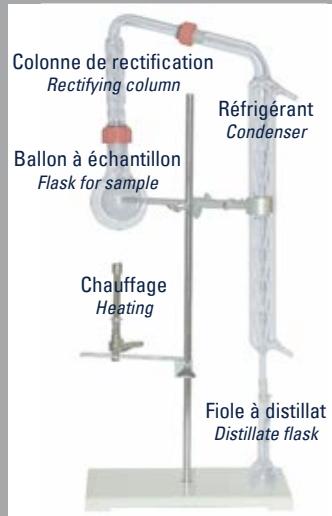
Alcohol is extracted by direct heating of sample in traditional glass stills or in automatic devices like Alco02. This method is recommended for beverages with high alcohol content (from 0 to 80% A.b.V.).

It is also possible to extract by steam the alcohol vapors (recommended method from 0 to 20% A.b.V., possible up to 40% A.b.V. with some conditions). In the same way, traditional models are available along with automatic devices DE EVO and DE2000. This same method is used for the extraction of volatile acids and sorbic acid.

A third range of instruments includes the traditional and automatic extractors like Vola 2000, exclusively designed for the extraction of volatile acids and sorbic acid by steam.

Determination method after distillation

- **Alcohol :** Density of collected distillate is linked to the A.b.V. It is measured by alcoholmetry, pycnometry, electronic densimeter or hydrostatic balance.
- **Volatile acidity :** Raw volatile acidity is determined by manual or automatic titration. The acidity of free and total SO₂ must be deducted to obtain the corrected volatile acidity.
- **Sorbic acid :** the concentration of sorbic acid in the distillate is determined by measuring the absorbance at 256 nm by means of a spectrophotometer, comparing it with the ones of a range of standard samples.

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Automatic distillers-extractors designed for the steam distillation of alcohol and the steam extraction of volatile or sorbic acids. All operations comply with European regulation (EEC) Nr.2676/90 and OIV standards described in the Compendium of International Methods of Analysis of Wines and Musts.

Measurement range (alcohol) : 0-40% AbV; from 40 to 80% AbV with 1/2 dilution.

Sample type : wines, spirits, fermented beverages, fruit juices and pastes, tomato sauces, vinegars, etc.

Characteristics :

- Automatic stop combined to the closing of the circulating cooling system
- Integrated steam generator (2800W)
- Bubbler with heating rings for an efficient steam extraction
- Easy drainage of the samples, even if not liquid
- Easy care and maintenance, easy access to the parts of the device (glass parts, steam generator)
- Short work time : 5 to 6 min by operation
- Power supply : 220V - 50 Hz
- Water supply : minimum pressure 2 bar, recommended temp. 18°C, minimum flow 12 l/min.
- Dimensions (HxWxD), weight : 93x46x40 cm - 30 kg

Supplied with glassware and chemicals needed for the distillation of alcohol and the extraction of volatile acids (except demineralized water).

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DE2000. Selection of the operating time with a timer.

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DE EVO. End of the distillation determined by a weighing system. Equipped with a cleaning gun

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ALCO 02, automatic direct distiller for spirits and beverages. It complies with the OIV recommendations in the Compendium of International Methods of Spirituous Beverages of Vitivinicultral Origin.

Equipped with an original preheating/heating system, it reduces significantly the time of distillation compared to classical laboratory stills and avoids alcohol losses.

Measuring range : 20-80% AbV.

Sample type : wines, liquors, spirits, etc.

Characteristics :

- Selection of the operating time with a timer
- Automatic stop combined to the closing of the circulating cooling system
- Heating power : 3200 W
- Constant display of the temperature of the refrigeration liquid
- Power supply : 220V - 50 Hz
- Water supply : minimum pressure 2 bar, recommended temp. 18°C, minimum flow 12 l/min.
- Dimensions (HxWxD), weight : 92x44x33 cm - 20 kg

Supplied with glassware and chemicals needed for the distillation of alcohol.

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The key points of a laboratory distillation :

- Volume to volume : the volume of collected distillate, after adjustment with distilled water, must be exactly the same as the original sample. The same volumetric flask is used to measure the original sample and collect the distillate. The distillate temperature must be the same ($\pm 2^\circ\text{C}$) as the sample temperature.
- The water running through the condenser must be cold ($14-16^\circ\text{C}$) to ensure an optimal condensation of the alcohol vapors, with a high flow (about 10-12 l/min).

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Mini Free Eau EVO. Designed to optimize, whatever the room temperature, the refrigeration of one distiller-extractor like DE EVO or direct distiller like Alco 02.

This compact cool bath is equipped with a thermostatic controller, a water tank and an immersed pump with stirrer. Thanks to its simple structure, it is very easy to install and fill.

Characteristics :

- Nominal cooling efficiency 1700W (Room temperature = 32°C, water at 15°C)
- Dimensions (HxLxP), weight : 55x40x75 cm, 51 kg (empty)
- Bath temperature : adjustable from -6°C to amb. temperature
- Power supply : 230V-50/60Hz

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Free Eau EVO. Adapted to optimize condensation with one or more DE or ALCO type distillers.

- High-power refrigerating unit (refrigeration power 3750W, absorbed power 2080W)
- Adjustable temperature : 5°C - Ambient temperature
- Circulating pump : 29 litres/min

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"Chercheur" alcoholmeter 0/100% AbV. Well designed instrument for a perfect vertical stability in the probe and a high accuracy for this kind of instrument. Division : 1% AbV. Length ≈ 300 mm.

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"Chercheur" alcoholmeter calibrated in 2 points, covering the 0/100 % AbV. range with 3 instruments. Division : 0,5 % AbV. Length ≈ 260 mm

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Small size alcoholmeter to measure 100 ml of distillate. Division : 1% AbV. Length ≈ 145 mm.

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Precision alcoholmeter calibrated in 3 points, covering the 0/100 % AbV range with 3 instruments. Division : 0,5 % AbV. Length ≈ 300 mm

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Class II precision alcoholmeter calibrated in 3 points, covering the 0/100 % AbV range with 5 instruments. Division : 0,2 % AbV.

	Longueur / Length = 310 mm	Longueur / Length = 290 mm
172000	0/20	172060 60/80
172020	20/40	172080 80/100
172040	40/60	172055 55/77

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Class II alcoholmeter calibrated in 3 points. These alcoholmeters comply with the rules defined in French laws. Division : 0,1% AbV. Calibrated at 20°C.

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Standard models, covering the 0/100 % AbV range in 10 instruments.

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The same, approved by a certificate of type examination, delivered with a certificate of initial verification to confirm the conformity of the instrument to the reference text, issued after comparison to a COFRAC instrument.

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Alcoholmeter kit from 173 series, with an inspected thermometer, in plastic case with protective foam.

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Rack for 6 alcoholmeters. Metal structure with Epoxy coating.

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Précision : $\pm 5 \times 10^{-5}$ g/cm³, 0,01% Vol.

Bench electronic density meter ALM-155. This densimeter has been especially developed for the enological application, taking advantage of technology and measuring cell of upper-end densimeters, but selecting the sole components for the specific needs of enological field. The Peltier thermostat fixed at 20°C and the specific software enables to offer this densimeter at an incomparable price on the market, with optimal value for money.

Sampling by built-in peristaltic pump (approximately 8 ml of sample), simple calibration (air/distilled water), easy cell cleaning, no maintenance needs, auto recording of the results.

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Précision : $\pm 5 \times 10^{-5}$ g/cm³, 0,01% Vol.

Bench electronic density meter DA-645. Measure of the resonant frequency oscillation. Touch sensitive digital screen. Equipped with Peltier thermostat. The accuracy provided by this instrument is perfectly adapted to oenology and beverages fields.

Data memory for 100 measurement methods, 4 modes of stability according to required accuracy, possible expression of density in concentration. Automatic calibration with air/water, or with standard solution. Minimal sample 1 to 2 ml, injected by a syringe (optional pump). Output : LAN for PC, USB to download data on flash drive. Output for printer (printer not included).

Measurement range : 0 to 3 g/cm³, alcohol from 0 à 100,00%.

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Jaulmes pycnometer. The density is calculated weighing a container, whose weight and volume at 20°C are known precisely, full of the liquid to be measured, taking into account the variations of air thrust on the pycnometer. The official tables give the alcoholic strength by volume according to the density.

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Iron test tube. The reaction created by the iron reagent causes a sediment, whose height in the graduated tube gives an idea of the iron content in wine.

198101 Kit complet à 1 tube / Complete kit with 1 tube

198102 Kit complet à 2 tubes / Complete kit with 2 tubes

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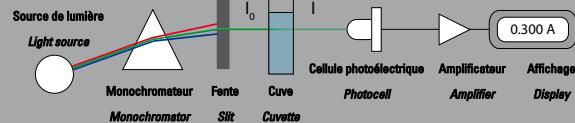
Spectrophotometry

Spectrophotometry is a quantitative and qualitative analytical method that consists of measuring the absorbance of a solution, i.e. its capacity to absorb light. The absorbance of a sample depends on the concentration of this sample and on the wavelength of the measurement.

Principle

When a light with intensity I_0 goes through a solution, a part of it is absorbed by one or more solute(s). The intensity I of the received light is lower than I_0 . The absorbance is defined by following relation : $A = \log(I_0/I)$. The obtained value does not have any unit. Beer-Lambert law defines the relation between wavelength, the solution absorbance, its concentration and the length of the light path.

Application

This principle is directly applied in oenology, to know the color intensity, which is the sum of absorbance values at various wavelengths (420, 520 and 620 nm), and to determine sorbic acid in a distillate obtained by steam extraction, after building a calibration curve with a range of standard samples.

Spectrophotometry is also used to determine easily a wide range of analytical parameters, useful for vinification and ageing, by means of enzymatic and colorimetric kits. The principle is to involve the compound to be determined in an enzymatic or colorimetric reaction that generates a chromophore chemical, which absorbs light at a given wavelength. Example : To know the malic acid content of a sample, reagents are added to induce the following enzymatic reaction : $L\text{-Malate} + \text{NAD}^+ \rightarrow \text{Oxaloacetate} + \text{NADH} + \text{H}^+$. Then the absorbance at 340 nm is measured, as NADH is chromophore at this wavelength.

We can offer a large range of spectrophotometers in order to meet user's needs :

- Analytical laboratories, with references 197200 and 197410, spectrophotometers with touch screen and extended features (spectrum analysis, concentration curves, etc.)
- Winemakers who need to control the most common analytical parameters with DS Oenophoto, a small and light photometer to perform one's own analysis at an affordable cost.

The main analytical parameters that can be measured with enzymatic and colorimetric kits are detailed p.36.

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DS Oenophoto. This photometer with several wave lengths enables fast and accurate analysis by the winemaker for common parameters such as residual sugars, yeast assimilable nitrogen (YAN), malic acid, etc. Completed with usual lab accessories and enzymatic or colorimetric kits, it gives an easy access to these analysis for all wineries.

Characteristics :

- Wavelength : 340 nm and 580 nm
- Absorbance : -0.2 à 3.0 A
- 1 position for macro or semi-micro cuvette with optical path of 10mm
- User-friendly app "DS Oenophoto" to download for free (control and data acquisition)
- Data exportation to an Excel file (csv file)
- Analysis up to 12 samples simultaneously
- Dimensions / weight : 11,5 x 6,5 x 5,5 cm / 230 g

Delivered with 3 micropipettes with disposable tips and a stand, 100 disposables cuvettes and a stand for cuvettes.

To be connected to a smartphone or tablet with Android (with Bluetooth connection), not supplied. Kits to order separately.

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Spectrophotometers 7200 (Visible) and 7205 (UV/Visible). They use diode array technology to produce exceptionally fast results. The colour touch screen interface provides fast and easy set up.

Characteristics :

- Modes : absorbance, concentration, kinetics, spectrum analysis in 50 points
- Small footprint and lightweight
- USB ports (communication to a PC and optional printer)
- Wide range of accessories available
- Equipped with a cuvette position 10x10 mm, and an external cuvette holder

Supplied with : internal memory 1 Go, 1 SD/USB drive 1Go, 100 disposable plastic cuvettes, user manual, power cable and software on CD-ROM with connection cable.

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Spectrophotometers 7410 (Visible) and 7415 (UV/Visible). They provide accurate and reliable results for various applications (industry, quality check...). The interface on a large 7" high definition colour touchscreen is fast and responsive.

Characteristics :

- Modes : absorbance, concentration, kinetics, spectrum analysis
- Android operating system
- Internal memory for data storage (10 Go)
- USB ports to store data on external support or to connect to printer (optional)
- Multilingual (french, english, german, spanish, italian)
- Extensive range of accessories available

Equipped with a single cuvette position 10x10 mm, in a wide chamber in which several accessories can be adapted.

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10 mm PS macro cuvettes (x100), volume 4,5 ml, disposable, for use between 320 and 1100 nm.

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10 mm PMMA macro cuvettes (x100), volume 4,5 ml, disposable, for use between 198 and 1100 nm.

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10 mm PS macro cuvettes (x100), volume 2,5 ml, disposable, for use between 320 and 1100 nm.

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10 mm PMMA macro cuvettes (x100), volume 2,5 ml, disposable, for use between 198 and 1100 nm.

- 5 197811 **Cuve en verre 10 mm.** Pour spectre visible.
10 mm glass cuvette, for visible spectrum.

- 6 197810 **Cuve en quartz 10 mm.** Pour spectre UV-visible.
10 mm quartz cuvette, for UV/visible spectrum.

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Other accessories for spectrophotometers : contact us.

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Thoma: Ruling with 16 big squares of 16 small squares each. Depth 0,1 mm. A big square = 4/1000 mm³.

Malassez: Ruling with 25 big rectangular fields with 20 small squares each. Depth 0,2 mm. A big rectangular field = 1/100 mm³.

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Growth media. Specifically conceived for analysis of wines and other fermented drinks. Pre-packages Petri dishes, available in 55 mm for samples on black membranes with grid, or 90 mm. Aerobic culture at 25°C in incubator (except Lactic Bacteria : anaerobic culture).

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Anaerobic system. For the incubation of one Petri dish in a bag.

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*Vinobrett™ par Veriflow. Thanks to PCR technology, the VINOBRETT™ system quantifies both active and VBNC *Brettanomyces bruxellensis* in less than 4 hours, with only a few minutes of manipulation. It enables an on-site proactive and affordable monitoring of *Brettanomyces*, even on young wines during fermentation.*

The system involves a specific thermocycler, the pre-aliquoted reagents in tubes, the proprietary buffers and the test cassettes.

- High sensitivity : from 10 cells/mL
- Matrix compatibility : juice, wine, lees, barrel rinsate, colony PCR, enrichment broth, samples from environment
- Simple preparation of sample
- Qualitative and quantitative assay
- Specificity : *Brettanomyces bruxellensis*

198400	Thermocycleur Veriflow <i>Veriflow thermocycler</i>
198401	Lecteur de cassette Veriflow <i>Veriflow cassette reader</i>
198450	Kit Vinobrett (24 tests) <i>Vinobrett Kit (24 tests)</i>
198451	Réactif PCR Veriflow (24 tests) <i>Veriflow Brett PCR reagent (24 tests)</i>
198452	Tubes Digest Veriflow (24 tests) <i>Veriflow digest tube (24 tests)</i>

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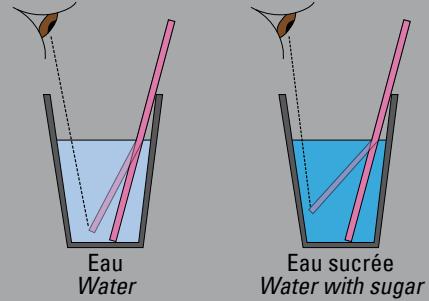
Refractometry

Principle

The refractive index is a characteristic of all liquid - or even pasty - media. The principle is : a light beam that goes from a medium to another changes direction, what determines a refractive angle, specific of the crossed medium. A homogeneous mixture of two components refracts light differently depending on the composition and concentration of this mixture.

The instrument used to determine the refractive angle, the refractometer, is composed of a prism where the sample is dropped and an optical or digital system for direct reading of the result. As the refractive index varies according to the sample temperature, a temperature correction, automatic in most cases, brings the readout to a value at the reference temperature (20°C).

Application

The refractometer is commonly used to determine the sugar content in simple media, such as fruit juices, musts, jam, etc. In this kind of sample, the other substances of dry matter have a very low influence on the refractive angle. The refractive index is converted so that the result is easier to read, generally in Brix degree, or in potential alcohol in oenology.

The refractive index is also used in many fields to control the quality of oils or the composition of antifreeze liquids for example.

Recommendations

The use of a refractometer is very simple and user-friendly, as the instrument is commonly equipped with a uniform sample distribution system on the prism.

How to take care of a refractometer :

- use preferably a plastic pipette to drop the sample on the prism ;
- ensure the device and the sample are at room temperature (difference between both below 5°C) ;
- check calibration on a frequent basis, by measuring distilled water or a standard solution ; adjust the 0 reading if possible ;
- clean the prism surface with water after use and dry with soft tissue ;
- avoid shocks or strong vibrations and store away from light and heat.

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Hand held refractometer Master TAL. This ergonomic refractometer is designed to easily and quickly measure the potential alcohol in the must with simplicity and reliability. Besides the automatic distribution of the sample, the instrument is equipped with a high quality of prism, an adjustable eyecup and an automatic temperature compensation from 10 to 30°C.

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RCT 0-22 refractometer. Water resistant and tough model to measure sugar content in grape, it is one of the essential item on the field. Equipped with an automatic temperature compensation, it gives reliable and accurate measures. Supplied with protective case. Range : 0/22% AbV Potential alcohol ±0,2.

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Hand-held refractometer RCT0-22 PA / 0-19 Baumé. With the same design as 196730 , it as a double-scale PA/Baumé dedicated to grape musts. Range : 0/22% AbV PA ±0,2, 0-19°Baumé ±0,2.

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Hand-held refractometer Master OE. Same design as Master TAL. It has a double-scale Oechsle/Brix to measure potential alcohol in must. Measuring range: 0/33% Brix - ±0,2%, 0/130° Oechsle - ±1°.

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Digital hand held refractometer PAL-Wine. Compact and water resistant, PAL-Wine is designed to measure the potential alcohol content in grape must. Calibration in accordance with the official method (E.C of 17/09/90). It is equipped with an automatic temperature compensation from 10 to 40°C, with a stainless steel sample stage and a wide liquid crystal display.

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Digital hand held refractometer PAL 84S, 86S, 87S. Same design as PAL-Wine, with specific scales.

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Digital refractometer WM7. It features 7 scales (AP, Brix, Oeschlé, KMW and Baumé.) and is developed to measure quickly and accurately sugar concentration in grape must. It is equipped with an automatic temperature compensation from 5 to 40°C, with a stainless steel sample stage and a wide liquid crystal display.

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Handheld refractometer, Master-alpha. Same design as Master-TAL. It enables the determination of sugar content in fruits, fruit juice and beverages. Measuring range : 0/32% Brix, ±0,2%

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RCT refractometers. Water resistant and tough models to measure sugar content in various sample types. Equipped with an automatic temperature compensation, they give reliable and accurate measures. Supplied with protective case.

	Désignation Designation	Gamme Range	Graduation Graduation
196729	RCT 0-18	0/18% Brix	0,2% Brix
196732	RCT 0-30	0/30% Brix	0,5% Brix
196734	RCT 5-54	5/54% Brix	0,5% Brix

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Digital pocket refractometers. Compact and water resistant, they measure the concentration of many different samples. Automatic temperature compensation from 10 to 75°C.

Désignation Designation	Gamme Range	Graduation Graduation
196711 PAL-1	0/53% Brix	0,2% Brix
196712 PAL-alpha	0/85% Brix	0,2% Brix

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RX5000i-Plus refractometer. This automatic digital refractometer is equipped with a sample cell on which the sample temperature is adjusted by Peltier effect. It displays quickly and precisely the refractive index, the Brix grade, or the concentration of many liquids. 100 user scales can be defined according to the samples to be analysed. Automatic recording of the 500 latest measurements. Temperature control range : 5 to 75°C Measuring range : 1,32420 to 1,58000 nD ±0,00002, 0,000 to 100,000 % Brix ± 0,001%

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pH METRY

pH is an essential tool for winemakers, as it is an indicator of acidity in musts and wine, and it helps to guide vinification. We offer a wide range of pH meters under several categories :

- *pH testers, which give an indication of pH but with low accuracy*
- *portable pH meters to get accurate results on the field*
- *bench pH meters, ideal for use at the laboratory, more ergonomic than portable version to be used on a table with stand for pH electrode*
- *pH meters and conductivity meters to measure conductivity and pH (conductivity meters are also available without pH parameter)*

Recommendations for use

When switching on the pH meter, calibrate the electrode with at least two buffer solutions (pH4 and pH7 for example). Buffer solutions must be thrown away after use. They have a short life after being in contact with air as pH changes quickly and calibration is no longer right. After each measurement, rinse the electrode with distilled water and dry it with a soft paper tissue before performing another analysis or before storage.

With a glass electrode, during a measurement, check that the ceramic junction is properly immersed (1), that the level of liquid in the stem is high enough (2), and that the stopper on the electrode is removed (3). Make sure this stopper is put back on the electrode after use.

Recommendations for storage

To avoid deposits and ensure a fast response, the bulb and the bulb junction with the stem must remain damp. If the pH meter is stored and after each use, clean the electrode and keep it in a storage solution (potassium chloride). Never store the electrode in distilled water. Store the electrode vertically, the bulb at the bottom. A frequent cleaning with a cleaning solution is recommended.

Used at room temperature and in good conditions, an electrode can be used for 1 to 3 years. An electrode not maintained properly or too old increases the response time, gives inaccurate results and requires more and more frequent calibrations.

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Checker electronic pH tester. Low cost electronic digital pH tester. Replaces pH litmus paper. Automatic calibration in 2 points. Range : pH 0 to 14. Accuracy (20°C) : ±0,2.

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GONDO pH tester. This waterproof pH tester is equipped with a digital display showing simultaneously pH and temperature. Automatic calibration in 2 or 3 points. Delivered in case with buffers pH4 and pH7, storage solution, and 4 1,5V batteries. Accuracy : ±0,05.

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PH7 Vio portable pH meter. Professional device with color display, with illumination and contrast. Only 4 keys allow the user to control all functions in an intuitive and easy way. Automatic calibration in 2 or 3 points.

This complete kit includes a pH electrode with integrated temperature sensor, colored pH buffer solutions, carrying case and accessories. Measurement range : pH 0 to 14, resolution pH : 0,1/0,01.

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Five Go portable pH meter. Ergonomic and intuitive, the FiveGo is a user-friendly quality instrument complying with usual requirements of accuracy. It is equipped with an automatic calibration in 3 points, an automatic temperature compensation and data memory for 200 readouts. Delivered with all accessories, ready to use, with a probe holder, a strap, a pH probe with temperature sensor and 2 buffer solution sachets (4.01 and 7.01).

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Bench pH meter pH50 Vio. This bench pH meter is both easy-to-use and accurate, with a very good value for money. It is equipped with an extra wide LCD with backlight showing simultaneously pH, temperature and measurement stability, and with an automatic temperature compensation. Delivered with combined pH electrode, temperature probe, electrode holder, buffer solutions sachets and electrolyte solution. Measurement range : 0,00 to 14,00 ; Accuracy: $\pm 0,01$.

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Bench pH meter Five Easy. Small foot print, equipped with a practical electrode holder; user-friendly : the FiveEasy is a precise and reliable device, with good value for money. Same technical features and electrode as FiveGo.

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COND7 Vio portable conductivity meter. Professional device with color display, with illumination and contrast. Only 4 keys allow the user to control all functions in an intuitive and easy way. Automatic calibration in 1 to 4 points.

This complete kit includes a probe with integrated temperature sensor, calibration solutions, carrying case and accessories. Measurement range : 0,1 μ S to 200 mS, accuracy : $\pm 2\%$ f.s.

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Five Go portable conductivity meter. Ergonomic and intuitive, it is equipped with an automatic calibration in 1 point, an automatic temperature compensation and data memory for 200 readouts. Delivered with all accessories, ready to use, with a probe holder, a strap, a probe with temperature sensor and 2 standard solution sachets.

Measurement range : 0,00 μ S/cm to 200 mS/cm ; Accuracy : $\pm 0,5\%$ of displayed value.

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710300 2

710200 3

Mallette du Seven Go Duo
Seven Go Duo carrying case

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PC5 pH and conductivity tester. Compact and waterproof, it is designed to measure pH, conductivity (EC) and Total Dissolved Solids (TDS) with automatic temperature compensation. Dual level display with 3-color backlight, showing measured value and temperature.

Measurement range : Cond 0-20 mS/cm, TDS : 0-10 ppt, pH -2 to 16, 0-50°C. Accuracy at 20°C : Cond $\pm 1\%$ f.s., pH $\pm 0,1, \pm 0,2^\circ\text{C}$.

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PC70 Vio portable pH meter-conductivity meter. Professional device with color display, with illumination and contrast. 6 keys

control all functions in an intuitive way. USB output (data transfer to PC or power supply). Automatic calibration pH in 1 or 3 points, conductivity 1 to 4 points. Data memory for 500 results. Complete kit in carrying case: 2 electrodes (pH and conductivity with integrated temperature sensor), colored pH buffer solutions, software for PC.

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Seven Go Duo portable pH-conductivity meter. User-friendly 2-way measuring instrument with a high reproducibility level thanks to automatic endpoint recognition, automatic calibration (pH in 3 points) and temperature compensation. Data memory : 99 readouts. Delivered with 2 electrodes with 2-meter cables, rubber case, calibration sachets, sampling flasks, clip for 2 electrodes and a compact carrying case.

Measurement range : 0,10 µS/cm-500mS/cm, pH 0-14, -5 to 105°C. Accuracy at 20°C : EC $\pm 0,5\%$, pH $\pm 0,01, \pm 0,1^\circ\text{C}$

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2.2.6. Solutions et accessoires / Solutions and accessories

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Electrode holder with flexible arm.

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Cleaning solution for pH electrode, 250 ml.

4 971500 Solution de conservation pour électrode pH 500ml. *Storage solution for pH electrode, 500 ml.*

5 971050 Solution électrolyte 30 ml / *Electrolyte solution 30 ml*

6 Solutions tampon / Buffer solutions

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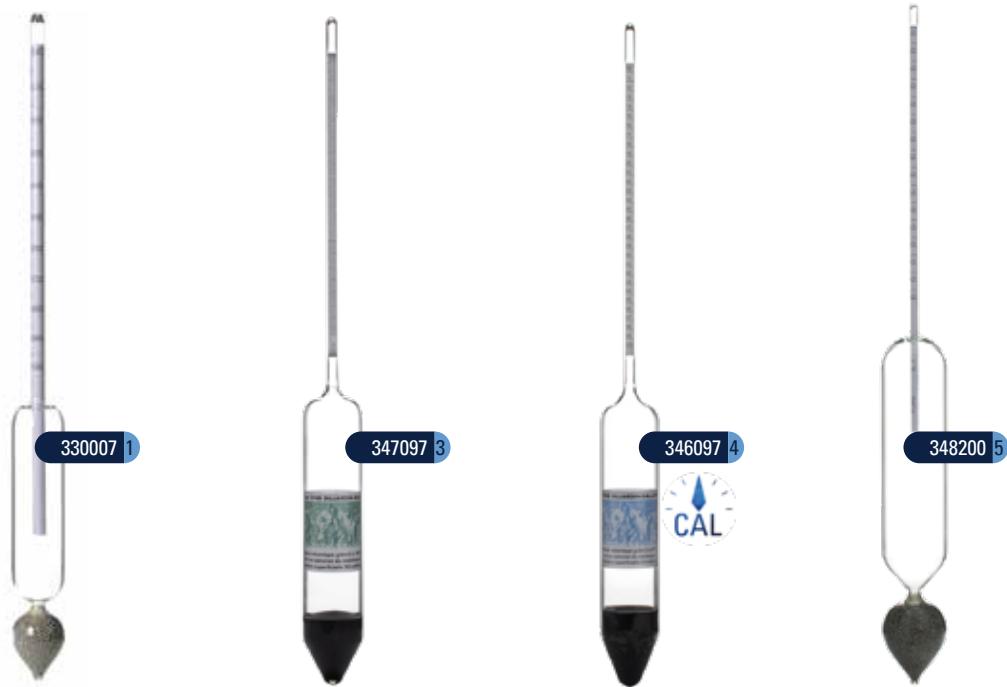
100 g range hydrometer, divided in 1 g/dm³. Length ≈ 280 mm

330007	700/800	330012	1200/1300
330008	800/900	330013	1300/1400
330009	900/1000	330014	1400/1500
330095	950/1050	330015	1500/1600
330010	1000/1100	330016	1600/1700
330105	1050/1150	330017	1700/1800
330011	1100/1200		

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Precision hydrometer : these hydrometers, covering 30 grams per instrument, comply with the rules defined in ISO norm. Division : 0,2 g/dm³. Calibrated at 20°C.

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The same, with an inspection certificate to confirm the conformity of the instrument to the reference text, issued after comparison to a COFRAC instrument.

Longueur / Length = 320 mm	Longueur / Length = 320 mm
346097	970/1000
346099	990/1020
346100	1000/1030
346103	1030/1060
346106	1060/1090
346109	1090/1120
346112	1120/1150

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Mustimeter with double scale (density 980-1130 g/dm³ and potential alcohol). According to EC rule Nr.2676/90. It allows to measure the sugar content in grape must according to its density and the potential alcohol content, to be obtained after complete fermentation.

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Guyot glucometer for grape musts, with triple scale (Baumé degree, potential alcohol and sugars in kg/hl). Length ≈ 300 mm

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Cider hydrometer. 1000/1100 hydrometer for apple must and ciders with double scale (density and potential alcohol).

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Beer hydrometer. Double scale (density 1000-1100 g/dm³, Plato degree) to check fermentation progress.

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Use of mustimeter : handling precautions

- The instrument must be perfectly clean and dry : clean with a solution of 20% sodium hydroxide, then rinse with distilled water and dry it with a clean tissue.
- Use it in a glass cylinder without graduation, straight (not conical). Do not use the instrument directly in the tipping trailer or tank.
- Make sure that the sample measured is representative. Remove rough particles to avoid wrong measurements (by settling, sieving, etc.).
- Immerse slowly the instrument in the sample, holding it at the top of the stem, giving an up and down movement, then release it with a rotary movement.

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Melli-mustimeter with triple scale (density, weight of honey in kg/hl and potential alcohol for honey wine)

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Hydrometers for various applications

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«Tirage» hydrometer 990/1005 g/dm³ divided in 0,1. Precise and sensitive model designed to be used for «tirage» wines. Calibrated at 15°C to be used in wine cellars. Length ≈ 325 mm.

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Extracto-enometer. Precision hydrometer to determine easily the dry extract in wines with hydrometric method (983-1003 g/dm³). Delivered with instructions and tables.

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Longueur ≈ 325 mm

Divided in 0,2 g/dm³, with an inspection certificate to confirm the conformity of the instrument to the reference text, issued after comparison to a COFRAC standard instrument. Length ≈ 325 mm

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Salinometer. Precision hydrometer to determine seawater density.

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Précision : ± 0,001 g/cm³ - 0,1°C

Portable electronic densimeter DA130N. This easy handling instrument, based on the measure of the resonance frequency, allows the measurement of density of many liquids, even viscous (wines, spirits, etc.). Automatic temperature compensation from 0 to 40°C. Output :IR port (Wire not included). Weight : 360 g

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Portable electronic densimeter Densito. With excellent ergonomics and an intuitive programme guided by a joystick, Densito offers fluid and fast process, with accuracy and repeatability. Automatic temperature compensation from 0 to 50°C. Output :USB port. Weight : 355 g

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The same, "Pro" version. Equipped with a RFID and bar-code to improve traceability of measurements.

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Red liquid filled laboratory thermometer

	Portée en °C Range in °C	Division Division
403009	-10/+60	1
403010	-10/+110	1
403007	-5/+35	1/2
403020	0/+35	1/2
403008	-5/+50	1/2
403021	0/+50	1/2
403022	-10/+50	1/2
403011	-10/+60	1/2
403023	0/+35	1/10

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PTFE protective sleeve. Application of a PTFE sleeve on a stem thermometer to protect marking from alcohols and to reduce the risks in case of breakage.

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Enclosed scale thermometer, red liquid, 0/+40°C, division 1/2.

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Red spirit filled thermometer with protective sleeve

	Portée en °C / Range in °C	Gaine / Sleeve
408000	-10/+60	Polypro
408003	-10/+120	Polypro
409000	-10/+120	Bois / Wood

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Use of an expansion thermometer

Before use, check that the liquid in the column is not separated due to rough handling in transport or incorrect storage. In case of separated liquid, it is necessary to reunite it by heating, cooling or centrifuging. Position the thermometer vertically immediately after this operation. Check that the thermometer is used in its measuring range. Store vertically.

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DSTemp. Digital Thermometer with high accuracy -20/+120°C with Pt1000 probe. Resolution : 0,01°C.

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Electronic thermometer for tank. K type thermocouple thermometer with LCD, -50/+300°C, ±1°C. HOLD function.

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Stainless steel probes for digital thermometer 411040 and 411010.

- 411039 30 cm, rigide / 30 cm, rigid
- 411101 150 cm, rigide / 150 cm, rigid
- 411102 210 cm, rigide / 210 cm, rigid
- 411103 150 cm, souple / 150 cm, flexible

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HQ1130 portable oximeter. Equipped with a LDO probe : no membranes to clean or replace, no electrolyte solution to refill, reduced calibration frequency, adapted to any kind of wine or beverage. The only maintenance is the replacement of the sensor cap only once a year. Automatic atmospheric pressure compensation. Data memory : 5000 readouts.

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Spare LDO lab probe for HQ30D oxymeter.

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Spare LDO field probe for HQ30D oxymeter.

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Bentotest, complete kit for the quick determination of proteic stability. It includes a 25 ml cylinder, reagents, a funnel and folded filters.

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Portable turbidimeter HI98703-02. Instrument with good value for money to determine turbidity during white wines vinification.

Nephelometric method ratio of scattered and transmitted light. 3 measuring modes (normal, average, and continuous). Tungsten lamp. Data memory : 200 records. RS232 and USB outputs.

Delivered in carrying case with cells, silicon oil, calibration set, 12V adapter and batteries.

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Calibration set (3 standard solutions + 1 test 0)

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Portable turbidimeter 2100Q. Reference instrument, user-friendly, to determine turbidity at all steps of vinification and before bottling. Ratio turbidimetric determination using a primary nephelometric light scatter signal (90°) and transmitted light scatter signal. The two-detector optical system compensates color in the sample, light fluctuation, and stray light.

Tungsten lamp. Data memory : 500 records.

Delivered in carrying case with cells, calibration kit STABLCAL, silicon oil. Measurement range : 0 to 1000 NTU (Accuracy $\pm 2\%$)

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Bench turbidimeter TL2300. Covering a wide range of turbidity, it performs accurate measurements based on a ratio of signals from 4 sensors. The intuitive user interface on large full color touch screen makes setup, calibration and measurements easier.

Tungsten lamp. USB Port.

Delivered with sample cells, StablCal and Gelex standard samples, silicone oil, and dust cover.

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The clogging index (CI) gives information about the wine filterability. Its determination requires a 2-bar pressure on 400 ml of wine filtered on a membrane, Ø25 mm, porosity 0,65 µm. It is based on the measurement of the time of filtration of successive parts of a wine volume at room temperature. A stainless tank is mounted on the filter holder, and it is completed with :

- at the top : a set mounted on the lid of the tank with a control manometer 0 to 4 bar, 1 draining valve and connector diam.6 for nitrogen or compressed air connection.

- at the bottom : 1 stainless steel valve ¼" with a removable DIN male connector to install the filter support disc.

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Dissolved CO₂

All wines, still and sparkling, contain CO₂ (carbonic gas). In particular, CO₂ plays an important part in still wine tasting. Vinification must take into account this parameter to assure that the CO₂ content during bottling is adjusted to the wine profile.

The determination of dissolved CO₂ in still wines is done easily with a Carbodoseur, giving quite accurate results if the measurement is performed in appropriate conditions, in particular at low temperature (sample below 10°C). Some more sophisticated devices enable to obtain fine-tuned values thanks to NDIR (Non Dispersive Infra Red) technology.

Pressure

In bottles of sparkling and semi-sparkling wines, excess pressure due to carbon dioxide can be measured with an aphrometer (pressure gauge). Their structure may vary depending on the closure type of the bottle, but most of them include a pressure gauge (manometer), a holding system and a piercing system.

The measurement must be done in the following way: the aphrometer is placed on the bottle and the fixing bracket positioned so that the instrument is stable on the bottle and there is airtightness between the inner space of the bottle and the aphrometer body. After piercing the crown cap or the cork with the aphrometer, the bottle must be strongly shaken until pressure is stable. Then the reading is done. In the case of permanent aphrometer, the device replaces the crown cap and is placed directly on the bottle ring. It allows to control the pressure increase all along the secondary fermentation.

The measurement must be done on bottles whose temperature has been stable for at least 24 hours. If the measurement temperature is different from 20°C, a correction must be applied, multiplying the measured pressure by the appropriate coefficient.

The aphrometer can also be used on bottles filled with still wines and closed with a cork, in order to check bottling quality. It is available for this use under a simple structure, without any holding system, and with a vacuum/pressure gauge. Bottles with screw cap require an aphrometer with fixing bracket.

To comply with OIV norm "Overpressure measurement of sparkling wines", the aphrometers must be regularly verified (at least on an annual basis) by means of a calibration bench.

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Carbodoseur : this is the simplest and most appropriate method to control dissolved CO₂ at the winehouse. The measure is done with a stoppered cylinder filled with the sample to analyze. Once shaken, the cylinder spills out a certain volume of the sample, proportionately with the CO₂ content. Average accuracy of 50 mg/l. Comes complete in a plastic case with calculation tables.

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CooRe Automatic CO₂ monitor. This device gives an accurate and immediate measurement of dissolved CO₂ concentration in wine. With user-friendly operations and fast measurements, it is appropriate for a lot of applications at the winery.

The built-in piercing system is adapted to several containers and does not require any specific strength. The protective door avoids spilling.

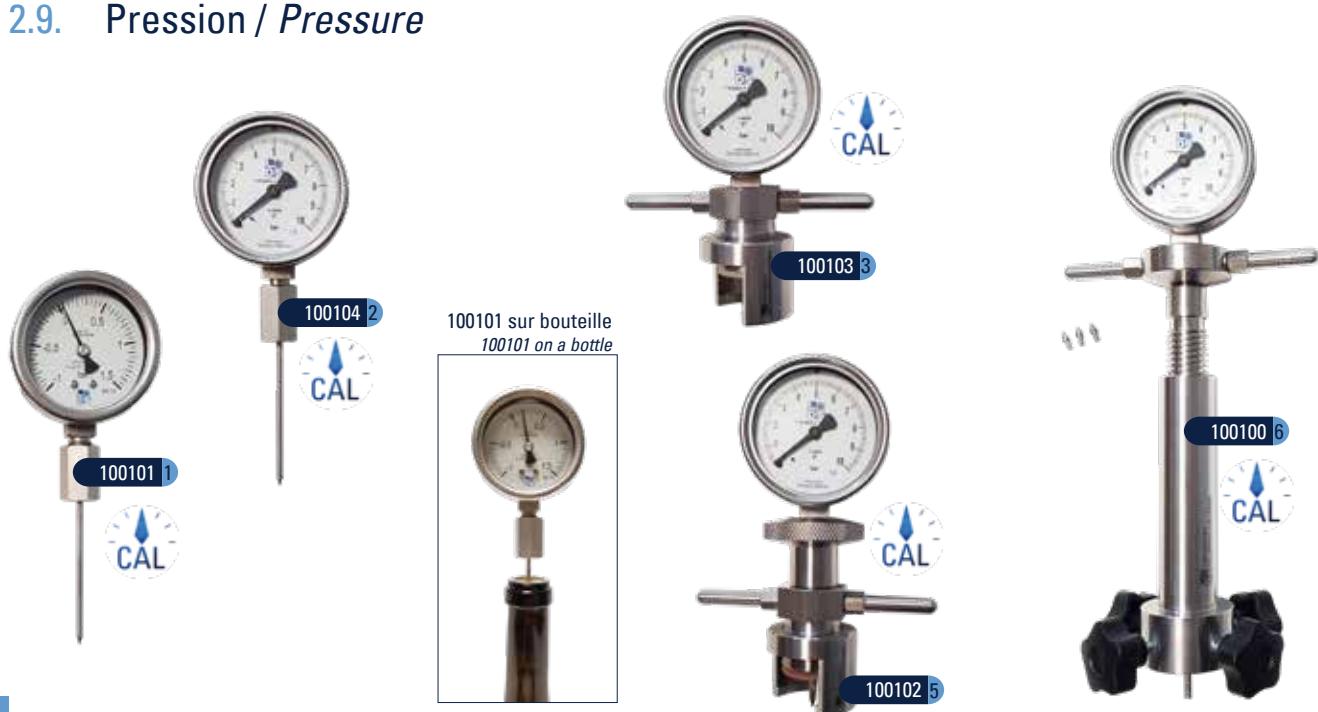
Characteristics :

- Control by touchscreen
- Analysis in 90 seconds, including piercing operation
- Measuring range : 0 to 24000 mg/L
- Temperature correction from 0 to 25°C
- Simple use, without specific training
- To be used with power supply or batteries, portable to carry it on site (12 kg)

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Simple aphrometer with manovacuometer -1/+1,5 bar to control the pressure and depression on bottles of still wine stoppered with cork.

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Simple aphrometer with manometer 0/10 bar to control the pressure on bottles stoppered with cork at the level of the neck.

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Permanent aphrometer with manometer 0/10 bar. Dedicated to permanent control of pressure increase during fermentation in bottle, it must be placed on the neck (Ø29mm) directly after bottling, without any cap or cork. Accuracy is guaranteed by a perfect airtightness of holding system.

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«Crown cap» aphrometer for sparkling wines. The 0/10-bar aphrometer shows the pressure instantaneously or permanently on bottle stoppered with crown cap (tinned iron and aluminum-stainless steel alloy). A gasket ensures airtightness after the needle goes through the crown cap.

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Cork» aphrometer for sparkling wines. Equipped with a 0-10 bar manometer, it allows to measure pressure instantaneously on bottles stoppered with cork. The detachable tip falls down inside the bottle after going through the cork to allow pressure measurement by the end of the needle.

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Options (integrated bidule, synthetic cork): contact us.

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Criotech. Quick and easy determination of tartaric stability based on electrochemical conductivity measurements under various conditions.

Two different methods are used : mini-contact and saturation temperature. Mini contact method determines KHT precipitation, expressed in μ Siemens. This value indicates the stability of the wine being tested. Saturation point : the analysis automatically provides the saturation temperature in °C of the tested sample.

Once the instrument has been prepared for an analysis, it proceeds automatically, requiring no further intervention by the operator until the last step (probe assembly cleaning).

Supplied with analytical software system to manage data : measurements readouts, in text or graphic mode (normal or differential), user status, single or multiple display mode : up to 10 analyses view simultaneously.

- Check control function allows easy and fast analyses.
- Touch screen display.
- Stepper motor guarantees repeatability of the measurements
- Silent ventilation system
- Automatic KHT dispensing
- USB interface to download data and upgrade the programme
- Standby function for energy saving

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Mini Logger USB is a single-use temperature recorder, in a waterproof pouch, identified by s/n, programmed to measure temperature every 11 minutes (maximal period of 60 days). It allows to control the temperature conditions during storage, shipment, etc. Comprehensive PDF report. LEDs on the key give information about overrunning the low (5 and 10°C) and high (20 and 25°C) temperature limits. Measurement range : -40 to +80°C; Accuracy : ±0,5°C Other settings : contact us.

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Hair tension hydrometer. Hygrometric scale 0 - 100%. Ø100 mm.

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Digital hygrometer-thermometer. Simultaneous display of hygrometry and temperature. MAX/MIN function. Measurement range : -10/+60°C (±1°C), 10 to 99% (±5%)

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Data logger for temperature and humidity KH70. Perfectly adapted to winehouses, this data logger is very simple to set up and read. Large LCD. 2 configurable setpoint alarms, with visual signal (LED). Magnetic mounting. Up to 50000 recording points. Lithium battery, approximately 500-day life (without display). Measurement range : 0 - 100 % RH / -20 to +70 °C

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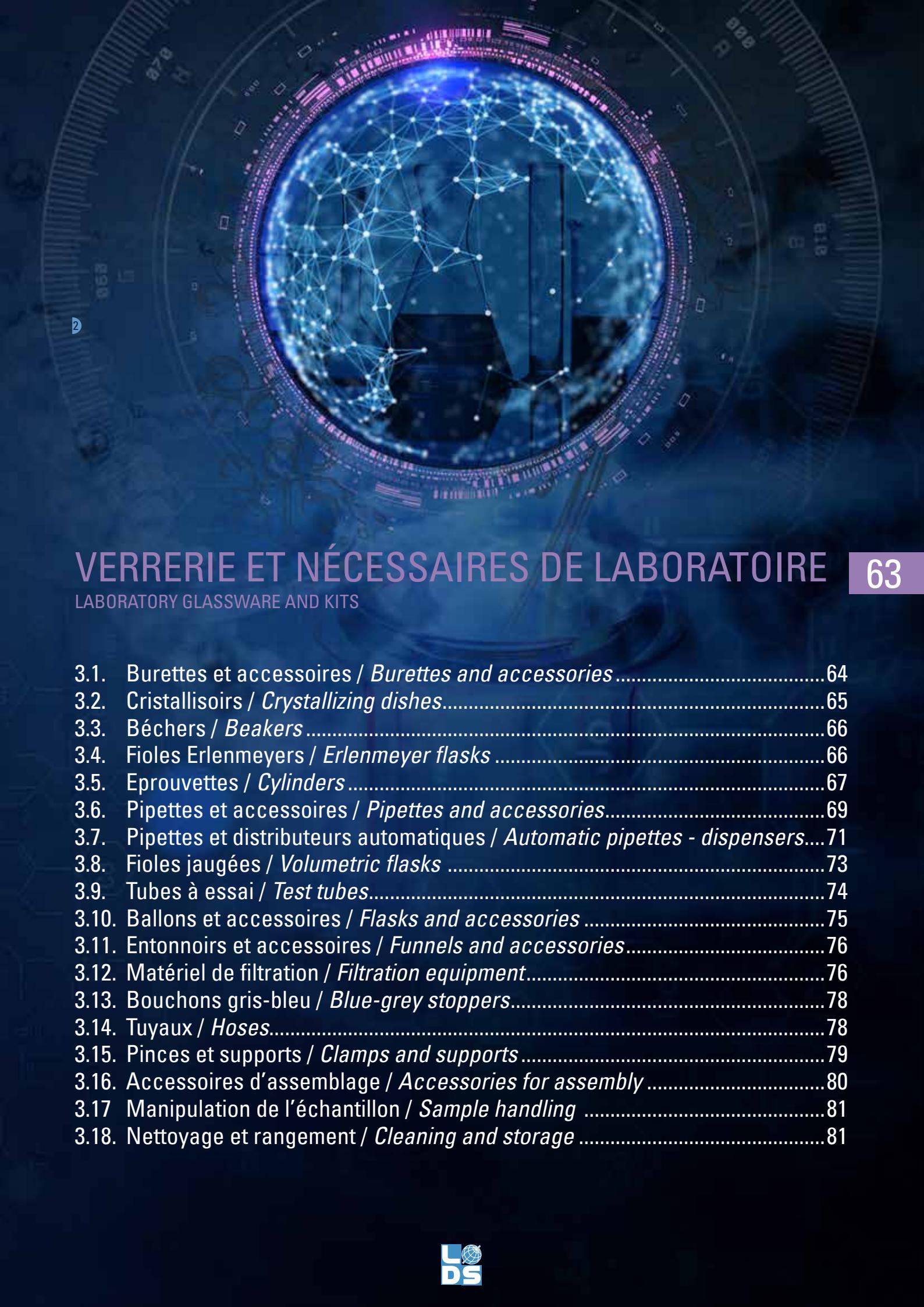
Anemometer with backlight LCD, waterproof, to be used at any time in any weather conditions. It became indispensable to respect safety requirements when spraying phytosanitary products.

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LABORATORY GLASSWARE AND KITS

3.1. Burettes et accessoires / Burettes and accessories

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Head for 24/29 ground joint flask for burette and bulb

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Schilling burette for a quick titration, automatic, mounted on a flexible PE flask and stand, equipped with pinch cock with titration screw for precise regulation of the flow.

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Burette with stopcock, engraved in mg/l of SO₂

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Burette with stopcock, engraved in acetimetric degrees

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- Pinces et supports / *Clamps and support*

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25 ml Brand bottle-top digital burette, piston-type, easy-to-use : rapid filling of the burette and precise titrations with the same hand wheels, 360° directional system. Easy and quick calibration. Quick air purge without air bubbles. Resolution 0,01 ml - Accuracy <0,1% Delivered without flask, with an injection tube, a telescopic filling tube, an assembly key, PP adapters and 2 AAA batteries.

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2,5 l amber glass flask for BRAND bottle-top burette

3.2. Cristallisoirs / *Crystallizing dishes*

3 Cristallisoir à bec en verre. DIN 12338.

Crystallizing dish with pouring spout, DIN 12338.

	Diamètre Diameter	Capacité Capacity	Hauteur Height
627040	40 mm	20 ml	25 mm
627050	50 mm	40 ml	30 mm
627060	60 mm	60 ml	35 mm
627070	70 mm	100 ml	40 mm
627080	80 mm	150 ml	45 mm
627090	95 mm	300 ml	55 mm
627100	115 mm	400 ml	65 mm
627150	140 mm	900 ml	75 mm
627170	190 mm	2000 ml	90 mm
627190	230 mm	3500 ml	100 mm

4 Cristallisoir sans bec en verre. DIN 12337.

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Crystallizing dish without pouring spout, DIN 12337. Diameter 70 mm, volume 100 ml, height 40 mm.

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Graduated glass beaker, with pouring spout, low form, ISO3819.

	Capacité Capacity	Diamètre Diameter	Hauteur Height
613050	50 ml	42 mm	60 mm
613100	100 ml	50 mm	70 mm
613150	150 ml	60 mm	80 mm
613250	250 ml	70 mm	95 mm
613400	400 ml	80 mm	110 mm
613600	600 ml	90 mm	125 mm
613800	800 ml	100 mm	135 mm
613901	1000 ml	105 mm	145 mm
613902	2000 ml	130 mm	185 mm

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Graduated PP beaker, with pouring spout, low form, blue marking, ISO7056.

	Capacité Capacity	Diamètre Diameter	Hauteur Height
613050P	50 ml	41 mm	60 mm
613100P	100 ml	51 mm	72 mm
613250P	250 ml	71 mm	95 mm
613500P	500 ml	87 mm	119 mm
613901P	1000 ml	109 mm	147 mm

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Graduated glass Erlenmeyer flask, narrow neck, ISO1773

	Capacité Capacity	Diamètre base/col Base/neck diameter	Hauteur Height
615100E	100 ml	64/22 mm	105 mm
615250E	250 ml	85/34 mm	140 mm
615500E	500 ml	105/34 mm	175 mm
615901E	1000 ml	131/42 mm	220 mm
615902E	2000 ml	166/50 mm	280 mm

4 Fiole Erlenmeyer col large en verre. Fiole graduée, DIN12385.

Graduated glass Erlenmeyer flask, wide neck, DIN12385

	Capacité Capacity	Diamètre base/col Base/neck diameter	Hauteur Height
615100L	100 ml	64/34 mm	105 mm
615250L	250 ml	85/50 mm	135 mm
615500L	500 ml	105/50 mm	170 mm
615901L	1000 ml	131/50 mm	280 mm

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Non graduated glass cylinders, tall form without pouring spout, especially designed for hydrometers and alcoholmeters.

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450 ml cylinder for mustimeters, with thermometer slot.

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325 ml calibrated cylinder for high degree alcoholmeters (173 and 174 series, above 50% AbV). Depth : 310 mm - int. Ø : 36 mm

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400 ml calibrated cylinder for low degree alcoholmeters (173 and 174 series, below 50% AbV). Depth : 330 mm - int. Ø : 39 mm

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450 ml calibrated cylinder for all alcoholmeters from 173 and 174 series.

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Graduated glass cylinder, class A, tall form, with pouring spout and hexagonal base, ISO4788.

	Capacité Capacity	Diamètre Diameter	Hauteur Height
623005	5 ml	12,9 mm	115 mm
623010	10 ml	15,5 mm	135 mm
623025	25 ml	21,3 mm	160 mm
623050	50 ml	26,0 mm	195 mm
623100	100 ml	31,3 mm	240 mm
623250	250 ml	40,7 mm	320 mm
623500	500 ml	53,2 mm	380 mm
623901	1000 ml	67,0 mm	465 mm
623902	2000 ml	85,0 mm	535 mm

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Graduated glass cylinder with ground joint glass stopper, class A, tall form, blue graduation and hexagonal base, ISO4788.

	Capacité Capacity	Diamètre Diameter	Hauteur Height	Rodage Ground joint
625100	100 ml	31,3 mm	240 mm	24/29
625250	250 ml	40,7 mm	320 mm	29/32
625500	500 ml	53,2 mm	380 mm	34/35
625901	1000 ml	67,0 mm	465 mm	45/40
625902	2000 ml	85,0 mm	535 mm	45/40

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Graduated translucent polypropylene (PP) cylinder, tall form, with moulded graduations, pentagonal base, and pouring spout, in accordance with ISO6706.

	Capacité Capacity	Diamètre Diameter	Hauteur Height
624050R	50 ml	25,5 mm	199 mm
624100R	100 ml	30,5 mm	249 mm
624250R	250 ml	41,5 mm	315 mm
624500R	500 ml	55,0 mm	361 mm
624901R	1000 ml	66,0 mm	439 mm
624902R	2000 ml	84,0 mm	531 mm

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Graduated PMP (polymethylpentene) cylinder, crystal clear, tall form, with blue graduations, pentagonal base, and pouring spout, in accordance with ISO6706.

	Capacité Capacity	Diamètre Diameter	Hauteur Height
626050	50 ml	26,0 mm	199 mm
626100	100 ml	31,0 mm	249 mm
626250	250 ml	41,5 mm	315 mm
626500	500 ml	55,0 mm	361 mm
626901	1000 ml	66,0 mm	438 mm
626902	2000 ml	84 mm	531 mm

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Volumetric glass pipette, Class AS, with safety bulb, ISO648.

1 trait	2 traits	Capacité	Hauteur
1 mark	2 marks	Capacity	Height
602001	603001	1 ml	310 mm
602002	603002	2 ml	330 mm
	603004	4 ml	350 mm
602005	603005	5 ml	400 mm
602010	603010	10 ml	440 mm
602015	603015	15 ml	455 mm
602020	603020	20 ml	510 mm
602025	603025	25 ml	520 mm
602050	603050	50 ml	540 mm

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Graduated glass pipette, 1/10, Class AS, with safety bulb, ISO835.

	Capacité	Hauteur
	Capacity	Height
604001	1 ml	350 mm
604002	2 ml	350 mm
604005	5 ml	360 mm
604010	10 ml	360 mm
604020	20 ml	450 mm
604025	25 ml	450 mm
604050	50 ml	720 mm

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Pasteur pipette. Disposable PE pipette with bulb and graduations. Capacity 3 ml, graduation 0,5 ml.

- 605103 Stérile, en emballage individuel
Sterile, individually wrapped
605203 Non stérile / non-sterile

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Universal pipette filler bulb with 3 valves.

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Semi-automatic pipette filler. New system with bulb, PTFE and silicon valves, allows to fill, adjust and dispense with only one finger. Adapted to all pipettes, from 0.1ml to 100ml.

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Pipette controller fitting any hand, working with glass or plastic pipettes. Adjustable aspiration/dispensing speed. Simple and fast operation. Rechargeable NiMH battery.

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Automatic pipette rinsing set. Made of high density polyethylene (PE-HD). Quick rinsing of pipettes, with continuous water refilling thanks to a siphon draining. It can be completed with a pipette basket.

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PE-HD pipette basket. To dip pipettes into the pipette washer. Height : 650 mm.

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PE-HD pipette jar. To pre-clean pipettes in cleaning agents. Height : 500 ml.

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Micro and macropipettes Acura are equipped with a large digital display for permanent volume visibility. Innovative concept of tip ejection. User friendly calibration system, light weight, ultra-soft, effortless activation, swift-set user calibration system, entirely autoclavable at 121°C. Available in fixed or variable volumes.

	Volume	Capacité / Capacity
605061	fixe / fixed	2 ml
605062	fixe / fixed	5 ml
605063	fixe / fixed	10 ml
605053	variable	20-200 µl
605050	variable	100-1000 µl
605051	variable	0,5-5 ml
605052	variable	1-10 ml

2 605080 Support pour 3 micro/macropipettes Acura
Work station for 3 micro/macropipettes.

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CLINIPET+. Fixed volume pipettes whose handle and mechanism have been especially designed for this series. The pipette features reduced pipetting forces and excellent accuracy and precision combined with highly durable, UV resistant and fully autoclavable body.

	Capacité / Capacity
605305	5 µl
605301	10 µl
605310	100 µl
605320	200 µl
605375	750 µl
605390	1000 µl

5 605300 Support pour 4 micropipettes Clinipet+
Work station for 4 Clinipet+ micropipettes.

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Calibrex dispensers with variable volume. With high tech materials and ergonomic design, this range is dedicated to precise dispensing of liquids from a large variety of bottles and flasks, with long lasting performance. Instant volume setting. Autoclavable at 121°C fully assembled. PFA glass coated plunger that enables trouble free distribution of salt solutions, weak and strong acids, as well as bases. Supplied without flask, with 3 PP bottle adapters

	Volume Volume	Divisions Division	Précision / Accuracy Vol.min. Vol.max.
600161	1 - 5 ml	0,1 ml	< 0,5% < 0,1%
600162	1 - 10 ml	0,25 ml	< 0,5% < 0,1%
600163	2,5 - 25 ml	0,5 ml	< 0,5% < 0,1%
600164	5 - 50 ml	1,0 ml	< 0,5% < 0,1%

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Volumetric glass flask, class A, with standard ground joint and PET stopper, ISO1042, DIN12664. Delivered with batch certificate.

	Capacité Capacity	Diamètre Diameter	Hauteur Height	Rodage RIN Stopper size
618050D	50 ml	50 mm	140 mm	12/21
618100D	100 ml	60 mm	170 mm	14/23
618200D	200 ml	75 mm	210 mm	14/23
618250D	250 ml	80 mm	220 mm	14/23
618500D	500 ml	100 mm	260 mm	19/26
618901D	1000 ml	125 mm	300 mm	24/29
618902D	2000 ml	160 mm	370 mm	29/32

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Volumetric glass flask, with one graduation mark, class A, especially adapted for laboratory stills.

	Capacité / Capacity	Hauteur / Height
619200	200 ml	180 mm
619250	250 ml	185 mm
619300	300 ml	200 mm

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Glass flask with 2 marks for sugar analysis, without stopper.

	Capacité / Capacity	Hauteur / Height
616050	50/55 ml	155 mm
616100	100/110 ml	180 mm
616200	200/220 ml	200 mm

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Very precise volumetric flask to check fill volumes. Factory calibration certificate not included, available upon order.

	Capacité Capacity	Graduation Graduation	Hauteur Height
617037	375 ml	365-385 ml	340 mm
617075	750 ml	680-780 ml	460 mm
617100	1000 ml	940-1040 ml	490 mm
617150	1500 ml	1440-1540 ml	520 mm

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Wooden protective case for calibrated fill flasks.

	Pour fiole For flask
617037C	617037 (375 ml)
617075C	617075 (750 ml)
617100C	617100 (1000 ml)
617150C	617150 (1500 ml)

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Automatic wooden tong. Length 250 mm. Grips a minimum diameter of 10 mm, maximum 35 mm.

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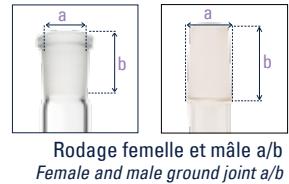
LABORATORY GLASSWARE AND KITS

3.10. Ballons et accessoires / Flasks and accessories

Rodage femelle et mâle a/b
Female and male ground joint a/b

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Round bottom glass flask, ground joint neck (Rodaviss).

	Capacité Capacity	Ø ext. corps Body outer Ø	Rodage Ground joint	Hauteur Height
610250S	250 ml	85 mm	29/32	135 mm
610500S	500 ml	105 mm	29/32	150 mm
610901S	1000 ml	131 mm	29/32	200 mm

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Other flasks with special neck : contact us.

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PP flask holder for round bottom flasks up to 10 litres. Ø160 mm. It supports hot flasks up to 120°C. Autoclavable.

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- Pince bois / Wooden tong
- Chauffe-ballons / Heating mantle

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Glass funnels with short stem, for filter paper, 60° angle, ISO4798.

	Capacité Capacity	Ø supérieur Upper Ø	Ø tige Stem Ø	Long.tige Stem length
363001	8 ml	35 mm	6 mm	35 mm
363002	15 ml	45 mm	6 mm	45 mm
363003	30 ml	55 mm	8 mm	55 mm
363004	115 ml	75 mm	8 mm	75 mm
363005	140 ml	85 mm	9 mm	85 mm
363006	190 ml	100 mm	10 mm	100 mm

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PP funnels with short stem, 60° angle, handle with eyelet.

	Capacité Capacity	Ø supérieur Upper Ø	Ø tige Stem Ø	Long.tige Stem length
363010	88 ml	75 mm	7 mm	55 mm
363011	222 ml	100 mm	8 mm	77 mm
363012	342 ml	120 mm	11 mm	90 mm
363013	817 ml	150 mm	15 mm	95 mm

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Double funnel holder. PP holder with epoxy base and stainless steel rod. For funnels with diameter 50 to 120 mm.

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Standard filter paper, smooth, medium filtration for general use in laboratory. Characteristics : 64 g/m², thickness 0,15mm, porosity 10-20 µm. Pack of 100 folded filters.

N°	Diamètre Diameter	Pour entonnoir réf. For funnel ref.
650300	00B	100 mm 363001, 363002
650310	0B	130 mm 363003
650311	1B	150 mm 363004, 363010
650312	2B	190 mm 363005
650313	3B	250 mm 363006, 363011, 363012
650314	4B	330 mm 363013
650315	5B	400 mm
650316	6B	450 mm
650317	7B	500 mm

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Quick filtration system. Complete kit : 10 ml glass syringe with Luerlock connector, a filter holder (Ø25 mm) and 100 filter disks.

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Filter disks (x100) for quick filtration system. Ø25 mm, por. 10-20 µm.

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Complete kit for quick degasification by suction of samples before analysis. It includes : a glass funnel filled with stainless steel wool, a 500 ml filter flask with stopper and a PP water-jet pump.

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PP water-jet pump, with sleeve nut R3/4", thread adapter R1/2", hose connection (olive 10/12mm diam.) and vacuum connection (olive 6/9mm diam.) removable with GL14 screwcap.

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Glass filtration unit. Complete kit with 300 ml funnel, sintered glass filter holder (porosity 10 µm), 1000 ml filter flask and aluminium clamp. Suction pump not included.

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3-position suction filtration unit. For membranes with diameter 47 or 50mm. 300 ml stainless steel funnels, stainless steel filter holder, aluminium forceps and PTFE valves. Suction pump not included.

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Filter membranes, made of mixed cellulose esters. Suitable for microbiology, colony counting, sample prefiltration and clarification. Diameter 47 mm, box of 100 units.

	Porosité Capacity	Couleur / type Colour / type
650420	0,45 µm	Noire quadrillée / Black with grid
650410	0,45 µm	Blanche unie / Plain white
650411	0,65 µm	Blanche unie / Plain white
650412	0,8 µm	Blanche unie / Plain white
650413	1,2 µm	Blanche unie / Plain white

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Other sizes, porosities, colors : contact us

6 Fiole à vide.

A tubulure et olive en verre.

Filter flask, with glass tube and olive.

Capacité Capacity	Ø corps Body Ø	Ø col Neck Ø	Hauteur Height
617250	250 ml	85 mm	35 mm
617500	500 ml	105 mm	35 mm
617901	1000 ml	135 mm	45 mm

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Conical grey-blue stoppers, with low taper ratio. They ensure a good airtightness and adhesion. Available without hole (P) or with hole (T).

N°	Diamètre <i>Diameter</i>	Hauteur <i>Height</i>
650102P/T	2 14/17 mm	21 mm
650103P/T	3 16/19 mm	23 mm
650104P/T	4 18/21 mm	25 mm
650105P/T	5 19/23 mm	28 mm
650106P/T	6 22/26 mm	30 mm
650107P/T	7 24/29 mm	32 mm
650108P/T	8 26/32 mm	34 mm
650109P/T	9 28/34 mm	36 mm
650110P/T	10 31/37 mm	41 mm
650111P/T	11 34/40 mm	43 mm
650112P/T	12 36/43 mm	45 mm
650113P/T	13 40/49 mm	49 mm
650114P/T	14 46/55 mm	51 mm
650115P/T	15 50/61 mm	56 mm
650116P/T	16 56/63 mm	61 mm
650117P/T	17 61/71 mm	65 mm
650118P	18 69/84 mm	66 mm
650119P	19 80/100 mm	67 mm
650120P	20 93/112 mm	88 mm
650121P	21 100/130 mm	85 mm

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Versilic hose, translucent, biologically inert. Complies with European Pharmacopeia chemical test requirements. Temperature of use : -50/+200°C. Possible peak at 220°C. Sold by the meter.

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Natural rubber hose. Possible sterilization in wet heat. Easy to fit on connectors thanks to its elasticity. Temperature of use : -35/+70°C. Sold by the meter.

4 617502 Tuyau à vide. Gomme souple rouge, diamètre 8/18 mm. Vendu au mètre. *Suction hose, made of red flexible rubber. Diameter 8/18 mm. Sold by the meter.*

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LABORATORY GLASSWARE AND KITS

3.15. Pinces et supports / Clamps and supports

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Stainless steel rod with threaded base, with nut and washer.

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Clamp with nut for tube and burette. Ni/Cr. Opening diameter 5 to 50 mm. Diameter of the rod < 18 mm.

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PP clamp with nut for 2 burettes. Diam. of the rod 10 to 14 mm.

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Complete kit for burette with one 170 x 110 mm base, a stainless steel rod 380 x Ø 8 mm and a clamp with nut.

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Clamp for stand without nut. Cork coated. Opening ≈ 40mm.

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Double nut for rods with maximum diameter 14 mm. Nickel.

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3-finger clamp with 2 tightening wheels. Chromium and zinc. Rubber covered jaws for a better grip.

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Open ring for stand. Ni/Cr. For funnels, flasks and iron gauze with ceramic gauze.

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- Accessoires de chauffage / Heating fittings

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Silicon PTFE sealing ring. For SVL sliding joint.

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Silicon sealing ring. For SVL sliding joint.

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Rodavis screw assembly. Complete kit.

650162	19/26 mm
650161	29/32 mm
650160	24/29 mm

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PP connector for plastic hose. To assemble with GL14 screwcap

650180	Droit / Right
650181	Courbé / Curved

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Laboratory spatula. Robust and economical spatula, made of glass fiber filled nylon. Autoclavable.

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3 594004 Pince inox pointe courbe 145mm.
Stainless steel forceps, curved end. Length 145mm.

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Glass weighing scoop, for standard use. With hollow stem to aid introduction of the powder into narrow neck vessels. The flat base ensures stability.

	Capacité <i>Volume</i>	Longeur <i>Length</i>	Matériau <i>Material</i>
594010	3 ml	70 mm	Verre / Glass
594011	6 ml	85 mm	Verre / Glass
594012	10 ml	100 mm	Verre / Glass

5 199720 Parafilm. Largeur 5 cm. Rouleau de 75 m.
Parafilm roll, width 5 cm, length 75 m.

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Standard drying rack, 55 positions. Made of PVC coated steel wire, with drip tray. Height 640 mm, width 360 mm.

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Hanging drying rack, 48 positions. Made of PVC coated steel wire. Height 500 mm, width 400 mm.

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Wall drying rack, 72 positions. Polystyrene. Equipped with 72 removable 100 mm pegs and a draining groove. Delivered with hanging hooks.

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Wall or hanging glass rack, chrome-plated. 5 rows for around 20 stemmed glasses in total.

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- Solution Glass Cleaner / *Glass cleaner solution*

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Magnetic mini-stirrer. Variable speed, from 100 to 1000 rpm. For small volumes (maximum 1 litre). ABS tray, 120x120 mm.

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Heating magnetic stirrer. White glass ceramic top with excellent chemical resistance, easy to clean. Cast aluminium body shaped for stability. Analogic adjustment of speed (maximum 2000 rpm) and temperature (maximum 450°C). Powerful stirring up to 15 litres.

- Dimensions of tray/heated area : 150 x 150 / 120 x 120 mm
- Fast heating and accurate temperature stabilization
- «HOT» warning light when the plate temperature is above 50°C, even if the device is off.

Fitting for retort rod (not included).

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Bench autoclave. Low footspace.

Permanent indication of the cycle steps.

Water level monitoring of the reservoir and sterilization chamber.

Water tank 4,2 litres.

Safety : automatic control of the water level in the tank before the cycle starts ; over-heating, safety valve and overpressure switch.

	Volume de la chambre <i>Chamber volume</i>	Puissance Power	Temp. Temp. <i>Temp.</i>	Opération Modes
606500	16 litres (Ø230 x 410 mm)	1,4 kW	121 ou/or 134°C	4, 15 ou/or 30 minutes
606501	16 litres (Ø230 x 410 mm)	1,9 kW	121°C	8 programmes pré-installés
606502	24 litres (Ø260 x 450 mm)	2,4 kW	ou/or 135°C	8 pre-registered programmes

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Vertical autoclave. Simple design, effortless opening and closing.

7 pre-registered programmes. Automatic standby operation.

Precise temperature control with additional pressure monitoring.

Electro-mechanical safety lock, safety valve, low-water interrupt, overpressure protection switch, overheating protection switch, circuit breaker, fault indicator (visual and audible).

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Compact and economical laboratory balance. Excellent value for money. Large LCD. «TARE» function. PVC weighing plate. 9V battery.

	Portée Weighing range	Lecture Readout	Plateau Weighing plate
601707	500 g	0,1 g	Ø 150 mm
601704	2200 g	1 g	Ø 150 mm
601708	5200 g	1 g	Ø 150 mm

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Precision balance. Stainless steel weighing plate. Delivered with power supply adapter, control weight and protective cover. RS232 output for printer or PC. Integrated programmes : parts counting, percentage determination.

	Portée Weighing range	Lecture Readout	Plateau Weighing plate
601700	200 g	0,01 g	Ø 105 mm
601702	400 g	0,1 g	Ø 105 mm
601705	2000 g	0,1 g	130 x 130 mm
601706	6000 g	1 g	150 x 170 mm

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Compact analytical balance. Modern all glass draft wind shield for optimal visibility of the weighing object. GLP/ISO record keeping of weight values with date, time and serial number. Higher weighing speed by switching off the last digit. Totalising of weights (manual/automatic). Possible homologation in legal metrology.

Alibi memory: paperless archiving of up to 100,000 weighing results. Internal memory for up to 999 weighing results, 1000 items or recipe ingredients, 100 container weights, 100 users.

Dispensing mode: by pressing one key you can set all relevant parameters for the dispensing process.

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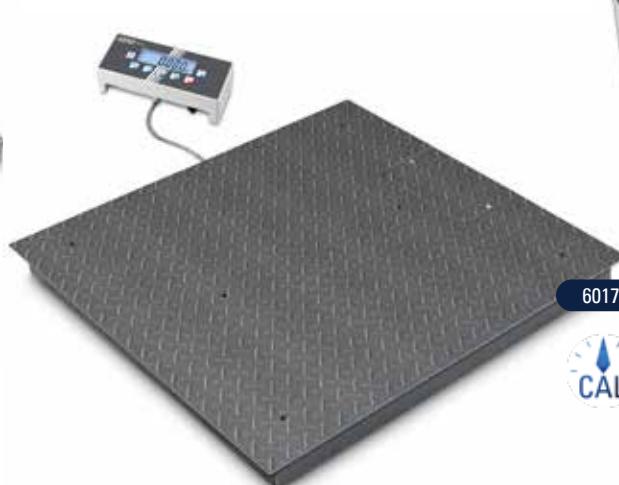
Scale with large platform. Simple and convenient 4-key operation. Wall mount for display device. Stainless steel platform, 55x55x6 cm. Power supply 230V or batteries.

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Floor scale. Weighing bridge made of anti-slip corrugated steel, 4 silicone-coated steel load cells. Easy levelling. Large display (25mm high figures). Universal power adapter. Optional ascending ramp or stable pit frame. Possible homologation in legal metrology.

	Portée Range	Lecture Readout	Charge min. Min. load	Plateau Weighing plate
601750	600 kg	0,2 kg	4 kg	120x150x11 cm
601751	1500 kg	0,5 kg	10 kg	120x100x11 cm

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Hanging scale. Perfect to weigh bags and buckets. Hold feature. Large display (12mm high figures). 3 AA 1,5V included batteries.

	Portée Weighing range	Lecture Readout
601760	50 kg	20 g
601762	100 kg	200 g

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DKD calibration (German calibration service). Valid internationally, it holds traceability to the national standard and determines the deviation from true, actual measure value.

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Compact ultrasonic bath. Capacity 0,75l. Digital control of timer : 90-180-280-380-480 sec. Bath opening (L x l x P) : 15 x 13 x 5 cm. Supplied with basket and lid.

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Ultrasonic bath. Indispensable at the laboratory to degas samples, to clean light pieces, etc. Rugged industrial transducers 40 kHz, a frequency that eliminates standing waves and creates an homogeneous cavitation in the tank.

Accessible control tower located behind and above the bath for safer operation.

Permanent operation/supply control, to maintain ultrasonic power, depending on the liquid level and temperature changes.

Drain valve with hose nipple and 1 m hose.

Plastic body, resistant to chemical products. Plastic tank cover included.

Mechanical timer, which can be set for up to 60 minutes and has a setting for continuous cleaning.

	Capacité Volume	Dimensions cuve (LxPxH) Tank dimensions (WxDxH)
609000	5,7 litres	29 x 15 x 15 cm
609001	9,5 litres	29 x 24 x 15 cm

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Additionally to the characteristics detailed above, this model is equipped with a 99-minute digital timer and continuous ultrasonic operation. High/low power control for more or less delicate applications.

	Capacité Volume	Dimensions cuve (LxPxH) Tank dimensions (WxDxH)
609005	5,7 litres	29 x 15 x 15 cm
609006	9,5 litres	29 x 24 x 15 cm

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PURA10 water bath. Equipped with a microprocessor maintaining a temperature stability of $\pm 0.15^\circ\text{C}$, this equipment offers many application possibilities. LED display and keyboard with splash-water protection. Easy-to-use. Lid not included.

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Transparent lift-up lid for PURA10 water bath, out of plexiglas and polycarbonate.

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SW22 shaking water bath. Equipped with a microprocessor for a $\pm 0.2^\circ\text{C}$ temperature control, this equipment is conceived for unsupervised operation : electronic timer, high/low temperature limits, automatic shut-down in case of low level. Splash-proof LED display and keyboard. Cover not included.

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Transparent lift-up Makrolon cover, for SW22 shaking water bath.

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Open Heating Bath Circulator C BT9. The circulator is mounted onto a bridge and can be easily removed. Transparent bath tank. Bright white display, easy to read. Precise temperature regulation, with stability $\pm 0,03^\circ\text{C}$. Very quiet and easy to use.

Working temperature : ambient to $+100^\circ\text{C}$.

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Refrigerated heating circulators to heat or cool external systems (refractometers, photometers, viscosimeters) and an internal bath tank made of stainless steel for small samples. High heating/cooling capacity to quickly reach the setpoint. Convenient cleaning of dust by removable venting grid. All models provide a drain port for conveniently draining the bath fluid.

- Bright white display temperature display, resolution 0.1 °C
- Lid included
- Temperature stability ±0,03°C
- Waste pump : 15 l/min, 0,35 bar.

Temp. de travail Capacité bain (LxPxH - volume)
Working temp. Bath capacity (WxDxH - volume)

608040	-20...+100°C	13 x 15 x 15 cm / 3...4 litres
608050	-40...+150°C	26 x 35 x 20 cm / 21...30 litres

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For all water baths, refrigerated and heating circulators with baths, many accessories are available (contact us) :

- Bath tanks to connect to circulators
- Tubes holder (different sizes)
- Adjustable platforms
- Bath fluids

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Dry heating block for 10 cuvettes. For the incubation of mixtures en cuvettes with optical path 10 mm before measuring the optical density. Temperature range : +25°C...+42°C (±0,1°C)

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Weight ring, for the stabilization of Erlenmeyer flask in an ultrasonic or thermostatic bath.

	Poids Weight	Ø intérieur Internal Ø	Pour fioles For flasks
606015	220 g	48 mm	125-500 ml
606016	450 g	56 mm	250-1000 ml
606017	680 g	61 mm	500-2000 ml
606018	910 g	74 mm	1 - 4 litres

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Ventilated bench centrifuge NEYA8. Very compact structure to optimize the space on laboratory benchtop. Quality raw materials, and stainless steel centrifugation bowl with easy access. Digital bright display and intuitive interface. Programmable centrifugation timer 30 sec. to 99:50 minutes. 3 acceleration and deceleration levels. Automatic locking system of the lid (motorized). Safety opening of the lid in case of absence of electric power. Induction motor, noiseless and maintenance free. Automatic rotor detection, check of presence of accessories and compatibility with maximum speed. Imbalance detection system with automatic stop.

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Ventilated bench centrifuge, NEYA10. With identical characteristics to NEYA8, it offers also the possibility to set 10 storable programs and to adjust speed in RCF. 9 acceleration and deceleration levels.

3 610111 Rotor libre SB4-175.

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Swing-out rotor SB4-175, including 4 B175 holders for NEYA8 or 10.

4 610115 Insert 4 tubes 15 ml for B175 holders (x4).

Tube insert for 4 15ml glass tubes for B175 holders (x4).

5 610121 Rotor angulaire A8-50.

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Fixed angle rotor A8-50. 8 holes for 50 ml tubes.

6 610123 Portoir et tube verre fond rond 50 ml.

Pour rotor A8-50.

Tube holder including 50 ml glass tube with round bottom. For A8-50 rotor.

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Complete range of rotors and tube holders : contact us.

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Glass alcohol lamp, 100 ml. With safe system, for wick Ø 4 mm.

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Flask holder, tripod, made of galvanized steel. Height approx. 21 cm.

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Iron wire gauze, made of galvanized iron wire mesh, with ceramic center.

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Gas burner with stopcock and pilot flame. Tube diameter 11 mm. For propane/butane gas.

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High flow burner Ø 40 mm. Ideal to heat big capacity water tanks (1000 ml).

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Gas burner holder with nut. It allows to adjust the height of the burner ref. 640123, fixing it on a rod with maximum diameter 12 mm.

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Metal-cased heating mantle. Plastic coated, chemical-resistant housing with built-in power-on and heating zone switch. Flexible glass yarn heating element. 2 heating zones (min/max). Built-in power regulator. Max. heating temperature 450°C. 1,5 m power supply (earthed) cable. Nominal voltage 230 V.

	Capacité Capacity	Puissance Heating power
606020	250 ml	150 W
606021	500 ml	200 W
606022	1000 ml	300 W
606023	2000 ml	500 W

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6-place heating mantle. Same structure as the above model, only one heating zone per position. Built-in power regulator with analogic adjustment at each position. Dimensions : 1150 x 150 x 290mm. Power supply : 230V.

	Capacité Capacity	Puissance Heating power
606050	250 ml	6 x 220 W
606051	500 ml	6 x 330 W
606052	1000 ml	6 x 495 W

3 606025 Noix pour tige verticale de diamètre maximum Ø12mm. Adaptable sur chauffe-ballons ref.60602X et 60605X.
Clamp for rod with diameter up to 12mm. Appropriate for heating mantles ref.60602X and 60605X.

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3-place heating mantle, or heating-mantles with other capacities, with magnetic stirrer: contact us.

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PH : 5.5 - 6.0

Dimensions - Poids : 50 x 15 x 45 cm - 10 kg

The BASIC PH4 water distillation unit is a continuous distillator. Easy to run, it is equipped with a safety thermostat. Production : 4 litres per hour thanks to a 3 kW chromium plated heater. The horizontal boiler is made of high quality borosilicate glass. CE certification and marking.

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Incubator CLN. Natural air convection. Stainless steel interior, plain door. Controlled by microprocessor. Digital display. Safety thermostat. Alarm in case of faulty temperature probe. Auto-diagnosis function. Control certificate at 37°C. RS232 output.

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Incubators with higher capacity, with refrigeration : contact us.

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Pocket magnifier with folding handle. Glass lens Ø40 mm. Chromium plated brass frame and handle. Magnifying factor: 4x

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Pocket lens made of organic glass. Black plastic case. Dimensions: 53x53 mm. Magnifying factor: 3,5x.

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Mini-microscope. This pocket instrument allows to visualize with precision the vine parasites thanks to an integrated lighting and to an appropriate magnifying factor (Magnifying factor : 30x).

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Stereomicroscope 20x/40x/80x, reflection/transmission type. Equipped with a 360° rotating head with 2x and 4x objectives. WF eyepieces 10x and 20x. LED illumination. Includes a dustcover, a pair of rubber eyecups, power supply adapter, glass plate and plastic plate (white/black).

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Researcher binocular microscope 40x-1000x. Transmission microscope adapted to school, study and laboratories. Nosepiece with DIN objectives 4x, 10x, 40x and 100x oil (achromatic). Eyepieces pair : WF 10x (Ø23mm). 20W halogen illumination with dimmer. Abbe condenser (N.A. 1.25).

Includes a dustcover, immersion oil and power cable. Trinocular model also available.

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Bioscience trinocular microscope 40x-1000x achromatic. Transmission microscope adapted to study and science. The ergonomic design facilitates its use. The trinocular head makes possible the simultaneous connection with the optional Mikrocam camera, for a permanent and simultaneous observation on screen. Nosepiece with DIN lenses 4x, 10x, 40x and 100x oil (achromatic). Eyepieces pair : WF 10x (Ø23mm). Coaxial mechanical desk with nonius scale. Coarse and precise focusing. 30W halogen illumination with dimmer, Kölher type. Abbe condenser (N.A. 1.25). Includes a dustcover, immersion oil and power cable.

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WF 20x eyepiece (Ø23mm), for Bioscience microscope.

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Mikrocam 1,3 MP 1280 x 1024. High quality of images. Possibility to make video sequences and single images. PC connection through USB 2.0 port. Professional software (supplied on a CD) with functions of image capture and measurement. Adjustable white balance and exposure control, manually or automatically. Supplied with 0.5x reducing lens, 23 mm adapter, and 2 rings 30 mm and 30.5 mm.

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Wide range of available accessories for microscopes : contact us.

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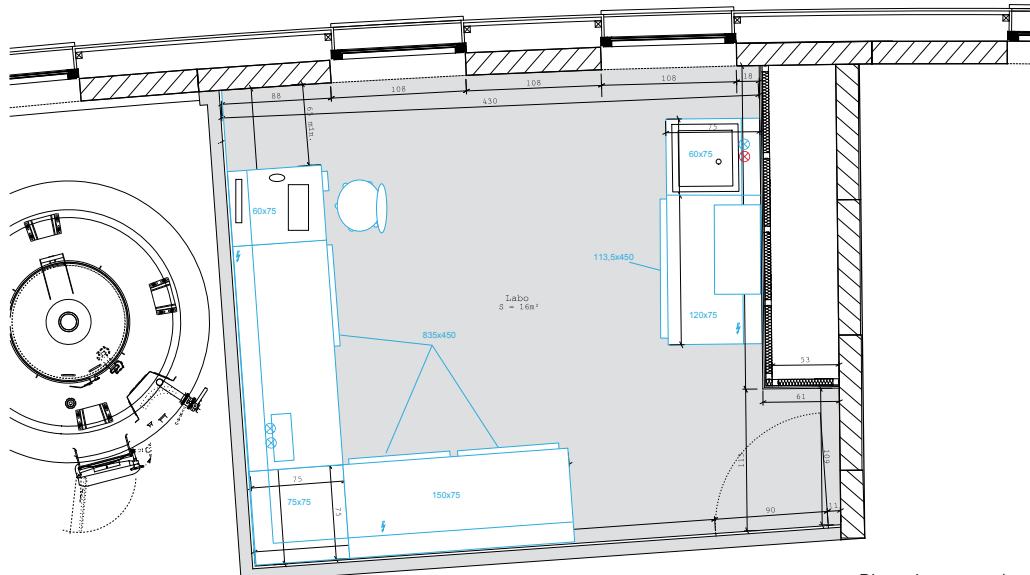
This furniture range offers a wide choice of dimensions, worktop materials, end pieces, fittings, etc. It has been conceived to be set up easily. The parts are delivered in spare parts but can be mounted upon request.

The elements hereafter are only a selection.

For an on-site or distance study, or to obtain information about the whole range (other coating/materials, other kind of furniture): contact us.

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16 m² laboratory approximately

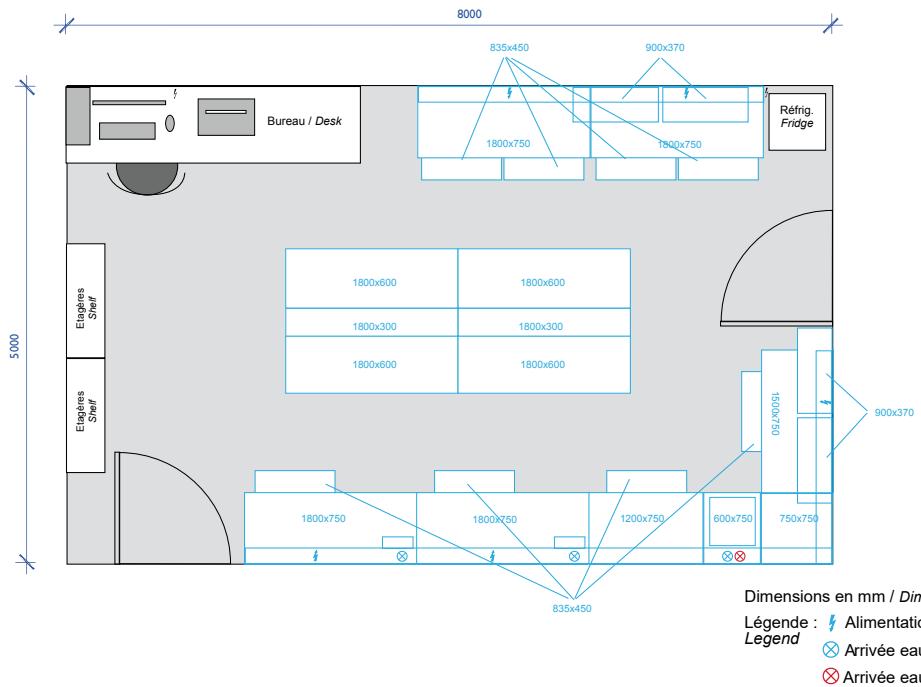
Suggested furniture (blue parts) :

- Dry benches with stratified work surface and top shelf on support, on metal leg
- Wet bench with stratified work surface and top shelf on support, on metal leg, equipped with a laboratory PP sink and epoxy-coated fittings
- Shelf under bench
- Stoneware sink unit with epoxy-coated fittings on furniture with top shelf on support
- Cabled 4 and 3-plug blocks
- Wall-mounted cupboard with solid doors
- Mobile cabinet with 4 drawers
- Mobile cabinet with 2 doors and a shelf
- Mobile cabinet with 2 doors and 1 drawer
- Laboratory chair on wheels
- Support for computer

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40 m² laboratory approximately

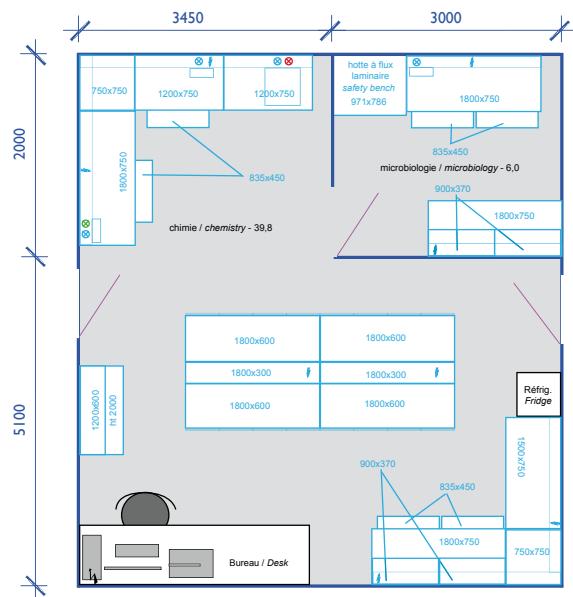
Suggested furniture (blue parts) :

- Dry benches with stratified work surface and top shelf on support, on metal leg
- Wet benches with stratified work surface and top shelf on support, on metal leg, equipped with a laboratory PP sink and epoxy-coated fittings
- Stoneware sink unit with epoxy-coated fittings on furniture with top shelf on support
- Central benches with technical box, stratified work surface
- Cabled 3-plug blocks
- Wall-mounted cupboard with solid doors
- Mobile cabinet with 4 drawers
- Mobile cabinet with 2 doors and 1 drawer

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45 m² laboratory approximately including :

- An area for chemical analysis with office section, 40 m²
- An area for microbiology, 6 m²

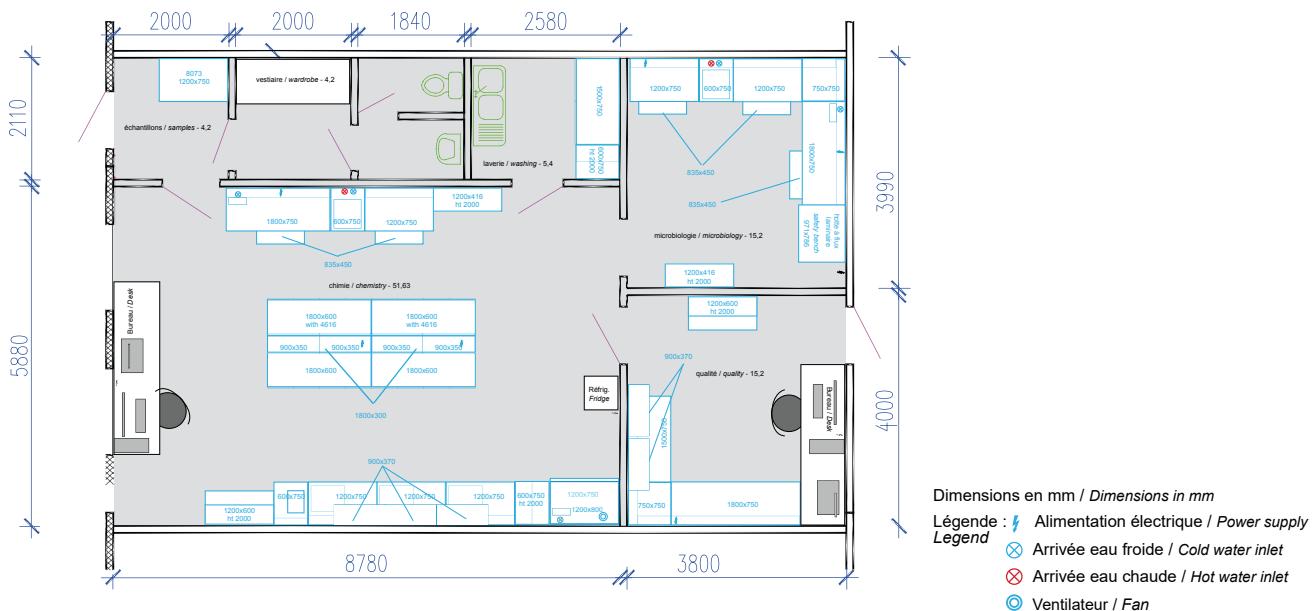
Suggested furniture (blue parts) :

- Dry benches with stratified work surface and top shelf on support, on metal leg
- Wet benches with stratified work surface and top shelf on support, on metal leg, equipped with a laboratory PP sink and epoxy-coated fittings
- Stoneware sink unit with epoxy-coated fittings on furniture with top shelf on support
- Central benches with technical box, stratified work surface
- Safety microbiological unit with laminar flow equipped with a UV lamp
- Cabled 3-plug blocks
- Wall-mounted cupboard with glass doors
- Mobile cabinet with 4 drawers
- Mobile cabinet with 2 doors and 1 drawer
- Built-in cabinet under bench with sliding doors
- 2-body cabinet (bottom with solid doors, top with glass doors)

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100 m² laboratory approximately including :

- An area for chemical analysis with office section, 50 m²
- An area for microbiology, 15 m²
- An area for quality control, 15 m²
- An area for washing, 6 m²
- An area for sample storage, 5m²

Suggested furniture (blue parts) :

- Dry benches with stratified work surface and top shelf on support, on metal leg
- Wet benches with stratified work surface and top shelf on support, on metal leg, equipped with a laboratory PP sink and epoxy-coated fittings
- Stoneware sink units with epoxy-coated fittings on furniture with top shelf on support
- Central benches with technical box, stratified work surface
- Laboratory fume hood equipped with a fan system
- Safety microbiological unit with laminar flow equipped with a UV lamp
- Anti-vibration table for analytical balance with top shelf on support
- Cabled 3-plug blocks
- Wall-mounted cupboard with glass doors
- Mobile cabinet with 4 drawers
- Mobile cabinet with 2 doors and 1 drawer
- Mobile cabinet with 1 drawer and 1 file drawer
- Cupboard with glass doors on both faces on technical box
- Built-in cabinet under bench with sliding doors
- 2-body cabinet (bottom with solid doors, top with glass doors)
- High cabinet with 2 vertical drawers
- Standard cabinet with 2 doors 4 shelves

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Multirisk safety cabinet, for the storage of non-flammable chemical products (acids, bases and solvents with low inflammability degree). Completely manufactured from 1,5 mm thickness steel sheet, cold pressed, acid resistant epoxy coated. One-piece structure with screw-jacks for level adjustment.

- Separation for "product category" in 3 compartments.
- Locking doors.
- Doors opening angle 110° to remove shelves without leaning them.
- Ring Ø125 mm for external connection or extraction system
- PP external hinges, corrosion resistant.
- Compartments equipped with basin form shelf, height adjustable, with welded corners.
- Safety signs complying with ISO 3864 norm.

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Cabinet with 3 solid doors, 3 compartments
Dimensions : 600 x 600 x 1950 mm

853003 Armoire 3 portes vitrées, 3 compartiments. Dimensions : 600 x 600 x 1950 mm
Cabinet with 3 glazed doors, 3 compartments
Dimensions : 600 x 600 x 1950 mm

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Accessories for multirisk safety cabinet.

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COMBISTORAGE acids/bases safety cabinet. Underbench version of multirisk safety cabinet. 2 compartments, 2 doors, 990x520x620 mm.

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Safety cabinet for flammable products. Firewall structure. Doors built with sleeves of various materials, to ensure a good resistance to fire. Resistant to high temperatures (800 °C).

- Automatic door closure in case the ambient temperature exceeds 50 °C ; Lock with key and door blockage ; Anti-sparking hinges.
- 30 mm thermo-expanding gasket (DIN 1402).
- Certified valves close the ventilation holes when the temperature exceeds 70 °C.
- Air extraction duct collar 100 mm diam. for a connection to outside or to a filtering aeration system.
- Grounding connection.
- Feet adjustment level.
- 3 epoxy coated shelves on stainless steel supports, height adjustable ; bottom spill pallet with draining grid.

853101 Armoire 1 porte, 3 étagères, bac de rétention et caillebotis. Dimensions : 595 x 600 x 1950 mm
Cabinet with 1 door, 3 shelves, retention container and spill pallet. Dimensions : 595 x 600 x 1950 mm

853102 Armoire 2 portes, 3 étagères, bac de rétention et caillebotis. Dimensions : 1200 x 600 x 1950 mm
Cabinet with 2 doors, 3 shelves, retention container and spill pallet. Dimensions : 1200 x 600 x 1950 mm

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Accessories for safety cabinet for flammable products.

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COMBISTORAGE safety cabinet for flammable products. Underbench version of safety cabinet for flammable products. 1 drawer, 990x520x620 mm.

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Accessory range, other extraction systems, other dimensions/ combinations: contact us.

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Dimensions externes (LxPxH) : 800 x 580 x 940 mm

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Vertical laminar flow cabinet. Conceived for microbiological and cell culture applications, it protects products or growth media (non-pathogen) in sterile environment.

Epoxy-coated steel frame. Perforated stainless steel (AISI 304) worktop, 2 panels, removable for cleaning. Temperate glass front window, possibility to lock in the fully open position.

Thermally protected motor fan, with manual regulation on control panel with "soft touch" membrane keyboard. Noise level: <58 dB(A). Absolute HEPA filter H14 class and pre-filters. Warning visual on display of exceeding the maximum period for the replacement of filters. Digital counter of working time.

Illumination work area > 700 lux (13W fluorescent white lamp)

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Classification : NFX15210 classe B.

Fume hood, EN "I" fire resistance class. Designed for chemistry laboratories, to protect operators from breathing harmful vapors of chemical products. Wood water/fireproof high-pressure laminate) panels and working area. Except for the metal lateral frame, acid-resistant epoxy coated. Plain sides. Back panel made of PP, removable for a total cleaning of the fume hood.

Low footprint on the bench. The side airfoil service panels are used for electric utilities on the left, and fluids on the right (water, gas, etc.). Mobile front sash. Made of Securit® glass 6 mm. Stop at 400 mm of the worktop. Break system on the sides.

Air flow optimization towards the centre of the fume hood to prevent vapors from going out.

Extraction system to be ordered separately.

Extraction outlet diam. 250 mm.

Total depth : 870 mm ; height 1600 mm.

Minimum height for installation : 2700 mm

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Fume hood with 1 front sash to be set up on existing bench.

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Floorstanding fume hood with 1 front sash. Metal support frame «C» shape, height 950 mm. Modular legs with epoxy-coated steel tubes. Adjustable screw-jack (minimum height 87 cm). Glass worktop. Optional back and side panels.

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Accessory range, extraction systems for fume hood, other worktop materials : contact us.

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Peroxyde test strips (x100). Enzymatic reagents. Results obtained by comparison with a colorimetric scale. Range 0-25 mg/l of H_2O_2 .

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Free and total chlorine portable photometer, wide range. LED light source, filter at 525 nm. Calibration control function. Intuitive use and measurements very easy to run. Delivered with 2 cuvettes and 3 1,5V AA batteries. Reagents not included.

Measurement range : 0,06 to 10,00 mg/l. Accuracy : $\pm 0,03 \text{ mg/l} \pm 3\%$ of reading ; excluding dilution error.

Free and total chlorine portable photometer, wide range. LED light source, filter at 525 nm. Calibration control function. Intuitive use and measurements very easy to run. Delivered with 2 cuvettes and 3 1,5V AA batteries. Reagents not included.

Measurement range : 0,06 to 10,00 mg/l. Accuracy : $\pm 0,03 \text{ mg/l} \pm 3\%$ of reading ; excluding dilution error.

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Reagent for 100 chlorine tests, wide range.

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Standard solution for free and total chlorine, wide range : 0,0 and 4,00 mg/l.

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Hardness (total) portable photometer. LED light source, filter at 466 nm. Calibration control function. Intuitive use and measurements very easy to run. Delivered with 2 cuvettes and 3 1,5V AA batteries. Reagents not included.

Measurement range :

Narrow range : 0 to 250 mg/L ; $\pm 5\text{mg/L} \pm 4\%$ of reading

Medium range : 200 to 500 mg/L ; $\pm 7\text{mg/L} \pm 3\%$ of reading

Wide range : 400 to 750 mg/L ; $\pm 10\text{ mg/L} \pm 2\%$ of reading

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Standard solution for total hardness.

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Lumitester Smart ATP-meter. This portable and light instrument gives practical information about the degree of hygiene of surfaces. The sample is taken by wiping a pad on a bounded surface, then this pad is inserted into a tube where it reacts with enzymes. This tube has to be inserted into the instrument to measure bioluminescence produced by this reaction.

Hygiene monitoring in three easy steps : Swab the sample, measure and analyse. Results within 10 seconds. Centralized multi-site data with cloud data management. Visualization and permanent control of data thanks to the application. Measurement in RLU – Relative Light Units. 2000 storage spaces

Delivered with 2 AA batteries, cleaning brush and USB cable.

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Lucipac Pen pads for Lumitester (x100). 5 packs of 20 tubes with enzymes and pads. To be kept cold (2-8°C) for one year.

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Sterile disposable patterns for ATPmeter, to define a 100 cm² area on a flat surface.

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Portable CO₂ detector. Light, ergonomic, high performance and with a long life battery (24h in continuous use), this detector allows people to work freely and stay in the winehouse and cellar without any risk.

Measurement by Nondispersive Infrared (NDIR) technology. Range : CO₂ 400-5000 ppm (± 30 ppm)

Definition of detection levels by the user.

Audible alarm in the event of danger, visible alarm (5 LED + LCD)

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Anti-scratch protective glasses. One lens piece. Can be worn over prescription glasses. Side air vent. Adjustment of temples in length and inclination

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Anti-scratch protective faceshield. Thermoformed anti-fog acetate faceshield. Adjustable elastic band, with soft foam anallergenic material at the front.

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Reusable earplugs. Pair linked by a cord. Soft and hypoallergenic material.

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Disposable particulate respirator with valve. FFP3 protection against particulates. Box of 20 pieces.

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Safety glasses for visitors. One lens piece. Can be worn over prescription glasses. Side air vent.

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HYGIEN AND SAFETY

6.3. Protection des personnes / Personal protection

199602 1

Raccord Easylock
Easylock connector

199611 2

Filtres non inclus
Filter not included

199616 3

Filtres non inclus
Filter not included

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Protection against pollution FFABE1P3 :

A1: protection against gas and organic vapors (boiling point > 65°C). Ex.: painting solvents, adhesives, pesticides.

B1: protection against gas and inorganic vapors (except carbon monoxide). Ex.: acetone, butane.

E1: protection against acids, sulphur dioxide (SO₂)

P3: protection against toxic dust and particulates, smoke, fogs (solid particles or aqueous aerosol). Ex.: copper, welding fumes, sodium hydroxide, calcium oxide, fibers, micro organisms, biochemical substances.

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Temporary half-mask (1-month use). FFABE1P3 protection, high performance. Integrated filters with flat shape for wider field of vision. Front piece made of thermoplastic material, hypoallergenic. Low resistance to respiration.

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Half-mask for Easylock filter holder. Front piece made of soft thermoplastic material, hypoallergenic. Equipped with a lace to let the mask hanging around the neck during breaks. Low maintenance.

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Complete mask with Easylock filter holder. Light conception. Frame with wide opening, easy to wear and remove. Easy maintenance. Panoramic lens with peripheral vision side panels.

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Preassembled filters : combination of a P3 dust filter and a A1B1E1K1 gas filter. For Easylock filter-holder. Box of 6 pieces.

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Latex, ambidextrous single-use gloves. Non sterile, powder-free, textured fingertips. Rolled cuffs minimise the risk of tearing. Suitable for food handling. Color: natural. Box of 100.

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Durable apron made from strong PVC. With nylon neck band. High resistance to oils, greases and certain chemicals.

Dimensions : 84 × 112 cm.

6.3.3. Premiers secours / First aid equipment

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Eye wash station, containing 2 solutions in flask to provide a hygienic and quick eye wash in case of contamination or irritation: «pH Neutral», a phosphate buffer solution for to neutralize acids and alkalis ; «Eye Wash», a sodium chloride solution to rinse dust and dirt. Ergonomic eye cup on the flask, for a hygienic and regular application.

	Forme Type	pH Neutral pH Neutral	Lave-œil Eye wash
199750	Station murale <i>Wall station</i>	200ml	500ml
199751	Coffret mural étanche <i>Waterproof wall case</i>	200ml	500ml
199752	Station murale <i>Wall station</i>	500 ml	1000 ml
199753	Coffret mural étanche <i>Waterproof wall case</i>	500 ml	1000 ml

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114

7.1.1. Tête-vin standards / Standard wine thieves

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Closed glass wine thief with rubber handle. Round end with a hole to be stoppered with the thumb. Equipped with a rubber handle, food grade.

Droit / Straight	Coudé / Curved	Longueur / Length
501025D	-	25 cm
501035D	501035C	35 cm
501045D	501045C	45 cm
501055D	501055C	55 cm
501065D	501065C	65 cm
501075D	501075C	75 cm
501095D	501095C	95 cm

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Personalized marking (one or more colors) : contact us.

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Open wine thief with rubber handle. Open end with a food grade removable rubber stopper with a hole to be stoppered with the thumb. Possibility to brush the inner surface with a bottle brush. Equipped with rubber handle, food grade.

Droit / Straight	Coudé / Curved	Longueur / Length
501125D	-	25 cm
501135D	501135C	35 cm
501145D	501145C	45 cm
501155D	501155C	55 cm
501165D	501165C	65 cm
501175D	501175C	75 cm
501195D	501195C	95 cm

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Straight wine thief with 4 holes, equipped with a rubber handle, food grade.

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Closed wine thief with glass handle(s). Round end with a hole to be stoppered with the thumb.

Droit / Straight	Coudé / Curved	Longeur / Length
504045	503045	45 cm
504055	503055	55 cm

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Glass tube for spirit sampling (topette).

507000	50 ml environ, diamètre 30mm 50 ml approximately, diameter 30 mm
507002	100 ml environ, diamètre 40mm 100 ml approximately, diameter 40 mm
507001	170 ml environ, diamètre 40mm 170 ml approximately, diameter 40mm

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Sample press. Adapted to fleshy fruit without stone. Equipped with a handle, aluminium structure, stainless steel cage and a drainpipe.

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Stainless steel ruler for barrel. Double graduation. With hanging ring. Length = 130 cm.

507100	En 1 partie / 1 part
507101	En 3 parties à visser / 3 parts to screw together

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Polypropylene jug, conical shape. Autoclavable at 121°C. Ergonomically designed handle.

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Aspirator bottle with spigot, made of high density polyethylene. Heavy duty material (2mm wall), conceived to be used for a long time. It includes a robust spigot to ensure a perfect leakproof performance. A few drops of vaseline shall be put on the thread of the spigot before assembly. 1 or 2 handles according to capacity.

	Capacité Volume	Diam.col Neck diam.	Diamètre Diameter	Hauteur Height
803001	5 l	46 mm	163 mm	335 mm
803002	10 l	55 mm	210 mm	414 mm
803003	25 l	55 mm	279 mm	525 mm
803004	50 l	55 mm	366 mm	603 mm

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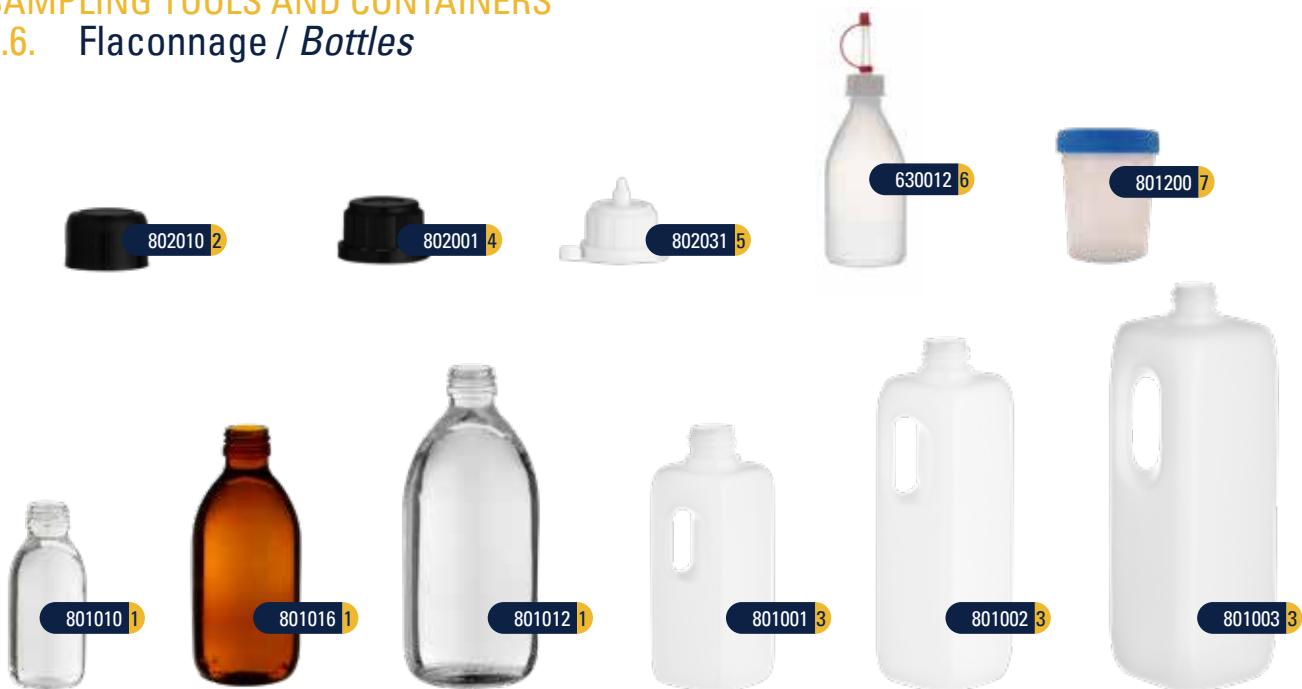
Tank with spigot. Made of high density polyethylene, food grade. Specifically designed to rationalize laboratory organization. 1 handle.

	Capacité Volume	Dimensions Dimensions
803011	10 l	380 x 300 x 160 mm
803012	20 l	470 x 370 x 180 mm

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7.6.1. Flacons en verre / Glass bottles

1 **Flacon type sirop en verre.** Bague PP28 (Ø ouverture 20 mm).

Glass flask, syrup type. PP28 ring (opening diameter 20 mm).

		Verre Glass	Capacité Volume	Diamètre Diameter	Hauteur Height
801013	Blanc / White flint	60 ml	38,6 mm	94,0 mm	
801010	Blanc / White flint	125 ml	49 mm	114,4 mm	
801011	Blanc / White flint	250 ml	61,7 mm	142,9 mm	
801012	Blanc / White flint	500 ml	76,8 mm	176,8 mm	
801014	Brun / Amber	60 ml	38,6 mm	94,0 mm	
801015	Brun / Amber	125 ml	49 mm	114,4 mm	
801016	Brun / Amber	250 ml	61,7 mm	142,9 mm	
801017	Brun / Amber	500 ml	76,8 mm	176,8 mm	

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Ribbed cap PP 28. For syrup bottle. Self seal.

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For other capacities or glass type : contact us.

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Invio 25 cap. Black, made of PE, with tamper-proof ring. For square bottle with handle.

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Invio 25 cap with dropper. White, made of PELD. For square bottle with handle.

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100ml dropping bottle, PE-LD, GL18 dropping insert and locking cap.

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PP cup. With blue screw cap. Capacity 150 ml.

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Wash bottle in polyethylene. Tube with dispensing tip which gives ultra-fine stream, removable to increase flow.

	Capacité Volume	Diamètre Diameter	Hauteur Height
601050	50 ml	35 mm	96 mm
601100	100 ml	43 mm	115 mm
601250	250 ml	60 mm	140 mm
601500	500 ml	75 mm	180 mm
601901	1000 ml	95 mm	220 mm

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Printed wash bottle, «Ethanol». Wide mouth with a valve that prevents liquid drips. Capacity 500 ml.

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Printed wash bottle, «Distilled water». Wide mouth. Capacity 500 ml.

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Our standard solutions and reagents are prepared and packed according to protocols described in the Quality Assurance Management System of the company.

The standard solutions are systematically controlled and adjusted before conditioning, using stabilized standard solutions.

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Acidimetric solution for total acidity.

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Acidimetric solution for volatile acidity.

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Acetimetric solution for acetic acid.

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Blue sodium hydroxide. With bromothymol blue. Concentration N/10. Short validity (1 month, 15 days after opening).

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Mixed solution for chromatography 250 ml. Short validity (1 month, 15 days after opening).

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931250 250 ml 931500 500 ml 931901 1000 ml

Soude (hydroxyde de sodium) / Sodium hydroxide.

Concentration 1,125N Concentration 2N Concentration 4N

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Colored reagent for Frantz Paul method, 500 ml.

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Milk of lime. Validity 2 years, 1 year after opening.

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Antifoam. Validity 1 year, 3-4 months after opening.

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Glass cleaner, ready-mixed alkaline cleaner for laboratory glassware.

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Titrivin offers a complete range of reference wines to calibrate apparatuses; and to assess and control methods of measurement or analysis. Complying with the ISO Guide 34, the batches are prepared by a COFRAC accredited laboratory for the activities of Reference Material Producer.

In order to guarantee the «wine matrix» effect, each reference wine is exclusively prepared by mixing wines and musts from various origins.

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The Titrivin AA series offers a range of ordered values of the 9 most common parameters to calibrate the response patterns of automatic equipments. Packaged in box of 20 glass vials of 10 ml.

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The Titrivin SW series offers a range of ordered values of the 9 most common parameters to calibrate the response patterns of automatic equipments. Increasing alcool content and decreasing sugars content, designing a matrix of fermenting musts or wines with residual sugars. Packaged in box of 20 glass vials of 10 ml.

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BTA Titrivin. Control standard for around 15 parameters. Used every 10 samples of wine analysed, it helps to control stability of measures on automatic analyzers. Box of 20 vials.

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BTB Titrivin. Control standard for manual analysis, with around 15 determined parameters. To be used periodically to validate the measurement and analysis methods run at the laboratory. Box of 6 bottles 250 ml.

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Ethykal, hydro-alcoholic solution with certified value of density. 250 ml flask, delivered with calibration certificate linked to a COFRAC certificate, giving the density and the calculated AbV. Approx. 12% AbV.

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Standard wine. Recommended for the calibration of electric ebulliometers. AbV determined by OIV official method. Certificate of analysis provided upon request.

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Malic acid, 3 standard samples for the determination of malic acid by chromatography on Whatman paper. 10 ml flask.

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Calibration solution for refractometer. 60 ml flask.

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Other concentrations up to 60% Brix : contact us.

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Practical alcoholometric guide, containing tables with :

- *Corrected alcohol content in % AbV according to the AbV reading and temperature. Range : -10 to +40°C (1/2°C) and 0 to 100% AbV (1/10 % AbV)*
- *Pure alcohol volume at 20°C in 100 dm³ of a hydro-alcoholic solution*
- *Content of methyl alcohol (% volume)*
- *Content of propyl alcohol*
- *Densities*
- *Calculations to increase/reduce the alcohol content of spirits.*

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Alcoholometric tables expressed in degrees. Alcohol content in % AbV according to the AbV reading and temperature. Range : 0 to +40°C (1°C) and 0 to 100% AbV (1% AbV).

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Gauging scale. Capacity of barrels, any shape.

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Inventory triple booklet. Capacity and volumes contained in containers. Extended range.

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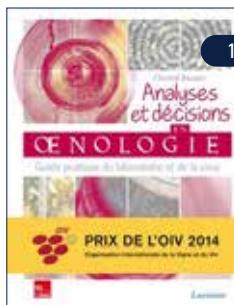
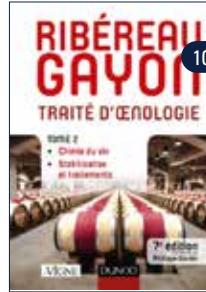
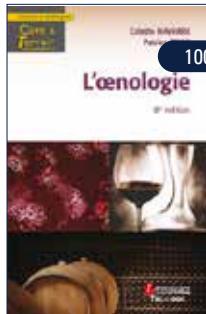
LIVRES

BOOKS

9.2. Livres techniques / Technical books

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Labox is a software with practical tools for operators in the field of alcoholic beverages (wines, ciders, beers, liquors, spirits, etc.). Thanks to a simple interface, it avoids the use of tables on paper and associated calculations.

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Software with alcoholmetric calculation : conversion of A.b.V. and volume as a function of temperature, recalculation of volume in pure alcohol.

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Alcoholmetry statement : configuration of cellars, positions and products for the inventory of pure alcohol

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Contact us for an online demonstration.

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Metrology is an essential part in product quality. It is a pillar in the same way as normalization and certification. As a matter of fact, accuracy of measuring instruments is fundamental to control manufacturing processes and assure the conformity of analyses.

Any measuring instrument is subjected to modifications due to wear and tear or clogging. Its metrological performances evolve, in a non-predictive way. Maintenance and regular control of measuring instruments allow to assure good results, to respect recommendation of ISO 9001:V2015 norm and to control procedures.

Laboratoires Dujardin-Salleron have set up the metrological department with the company E.R.IN (Study and Research in Industry), with the purpose of calibration, control, verification and maintenance of instruments. The provided services comply with norms in force and lead to issue a certificate that guarantees traceability to international system of units.

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REFRACTOMETRY : FROM 0 TO 72,8% BRIX

Type : handheld refractometer, optical or digital

Recommended frequency : on an annual basis, or after a long period without use

Self-control : distilled water or standard solutions (cf.p.125)

Metrological services : calibration and adjustment in 2 points

Method :

- use of a standard block, glass sample measured with the method of grazing incidence on a very accurate refractometer (Abbe refractometer)

- use of calibration solutions, manufactured and controlled internally following a normative procedure

THERMOMETRY : -20 to +150°C

Type : expansion (stem or enclosed type) or digital

Recommended frequency : on an annual basis for electronic thermometers, and according to internal norms of the user institution for expansion thermometers

Other occasions : if the thermometer has not been calibrated for several years and shows a drift

Metrological services :

- control in 1 point
- verification in 1 point (+0°C or initial value)
- calibration in 1 point (+0°C or initial value)
- verification in 3 points (+0°C or initial value)
- calibration in 3 points (+0°C or initial value)

Method : calibration bench consisting of a thermostatic bath and a traceable chain Pt100 4 wires

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Metrological services : verification and calibration

Method : calibration in 3 points (ambient air, zero, and air saturated with reference gas)

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HYDROMETERS – ALCOHOLMETERS

Type : glass hydrometers and alcoholmeters

Recommended frequency : according to internal norms of the user institution

Metrological services :

- simplified hydrostatic method, in 3 points, at 15°C or 20°C, uncertainty 0,04 g.dm³ + 0,3 division
- complete hydrostatic method, in 3 points, at 15°C or 20°C, uncertainty 0,02 g.dm³ + 0,1 division if length between 2 marks > 1 mm

Method : Cuckow calibration bench

Calibration of glass hydrometers requires to know the manufacturing or use characteristics of the instrument : the surface tension of the liquid that the hydrometer is supposed to measure, and the reading mode of the meniscus.

PRESSURE : -1 TO +15 BAR

Type : Bourdon tube pressure gauge

Recommended frequency : on an annual basis

Metrological service : calibration in 5 or 10 points, depending on the class and the diameter of the pressure gauge

Method : calibration bench consisting of a pneumatic pressure generator and a digital manometer connected to national standards

MICROPIPETTES : FROM 200 µL TO 10 ML

Type : piston-operated volumetric apparatus

Recommended frequency : according to internal norms of the user institution

Metrological service : verification in 1 point

Method : gravimetry according to ISO 8655 norm

VOLUMETRY : FROM 200 µL TO 10 L

Type : laboratory glassware, graduated or gauged

Recommended frequency : according to internal norms of the user institution

Metrological services :

- control in 1 point with mercury or water
- verification at the mark with mercury or water
- calibration at the mark with mercury or water
- verification in 3 points with mercury or water
- calibration at the mark with mercury or water

Methods :

- graduated cylinders according to NF EN ISO 4788
- 1-mark volumetric flask according to NF EN ISO 1042
- burettes according to NF EN ISO 385
- laboratory glassware according to NF EN ISO 384
- volumetric instruments according to NF EN ISO 4787

OVENS : 0 TO 1600°C

Type : any kind of oven

Recommended frequency : on an annual basis, or after a maintenance or repair operation

Metrological service : characterization of the oven

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ALPHABETICAL INDEX

A

Acetaldehyde 33
 Acetic acid
 Kit for titration (vinegars) 12
 Spectrophotometry 33
 ALCO 02 26
 Alcohol
 Alcoholmetry 28
 Distillation 23
 Ebulliometry 21
 Pycnometry 30
 Alcohol lamp 92
 Alcoholometers
 "chercheur" 28
 class II verified in 3 points 28,29
 in case 29
 precision verified in 3 points 28
 small model 28
 with certificate 29
 Alcoholometric tables in degree 128
 Anaerobic culture 38
 Anemometers 62
 Anthocyanins 34
 Aphrometers 60
 Apron 112
 Ascorbic acid 33
 Aspiration/Oxydation 17
 ATP-meter 109
 Autoclaves 85

B

Balances 86
 Beakers 66
 Benches 98
 Bentotest 56
 Books 129
 Botrytest 31
 Bottles 117
 Box of 12 wicks 21
 Brushes 81
 Buffer solutions 48
 Bulb for auto-zeroing burette 64
 Burettes
 auto-zeroing 7,64
 digital 65
 graduated in acetimetric degrees 65
 graduated in mg/l of SO₂ 64
 Mohr 7,64
 Schilling 64

C

Calibration 132
 Caps (screw assembly) 80
 Carbodoseur 59
 Centrifuges 91
 Ceramic gauze 92
 Chemical products 123
 Chromatography on paper 12

Cider hydrometer 50
 Citric acid 33
 Cleaning solution 48
 Clogging index 57
 CO₂ portable detector 110
 Conductimeters 46
 Conductivity standards 48
 Cooling bath with circulation 27
 Copper 33
 Criochek 61
 Crystallizing dishes 65
 Cuves for spectrophotometry 35
 Cylinders
 calibrated 67
 graduated 68
 without graduation 67
 with thermometer slot 67

D

DE 2000 25
 DE EVO 25
 Degassing kit 77
 Dissolved CO₂ 59
 Distillers-Extractors
 ALCO 02 26
 DE2000 25
 DE EVO 25
 Mothe style 23
 Dispensers 72
 Dropping pipette 69
 Drying racks 81
 DS Cenophoto 33
 Dynamometers 87

E

Earplugs 110
 Ebulliolog 22
 Ebulliometers
 Traditional ebulliometer 21
 Ebulliometer 21
 Ebulliotronic 22
 Electrodes
 pH 17, 48
 platinum 15,17
 Electrolyte solution 48
 Electronic density meters 30
 Enzymatic and colorimetric kits 36
 Erlenmeyer 66
 Ethycal 125
 Extraction apparatus for V.A.
 DE2000 25
 DE EVO 25
 Duclaux-Gayon style 9
 Mothe style 9,23
 Vola 2000 10

Extracto-œnometers 51
 Eye wash station 112

F

Fehling method 19
 Fermentest 19
 Ferrometer 31
 Filter membranes 77
 Filter stand 77
 Flask holders 75,92
 Flasks
 Erlenmeyer 66
 fill flask 74
 filter flask 77
 flat/round bottom 75
 volumetric 73

Flow detector 78
 Flowmeter for Frantz Paul 17
 Folded filters 76
 Forceps 74,79,81
 Frantz Paul apparatus 17,18
 Free Eau Evo 27
 Fume hood 105
 Funnels 76
 Funnel holder 76
 Furniture 98

G

Gas burners 92
 Gauging scale 128
 Glass micropipette 69
 Gloves 112
 Gluconic acid 33
 Glucose 33
 Glycerol 33
 Growing media 37
 Guyot glucometer 50

H

Heating mantles 93
 Hoses 78

Hydrometers
 15 grams range 49
 30 grams range 49
 60 grams range 49
 100 grams range 49
 Baumé 51
 extracto-œnometer 51
 for beers 50
 for ciders 50
 "tirage" 51
 with certificate 49

Hygrometers 62

I-J

Incubators 94
 Inventory triple booklet 128
 Iron
 Kit for determination 31
 Spectrophotometry 33

ALPHABETICAL INDEX

L

- Labox 131
- Laccase activity 31
- Lactic acid
 - Chromatography on Whatman paper 12
 - Spectrophotometry 33
- Laminar flow cabinet 104

M

- Macropipettes 71
- Magnetic stirrers 84
- Magnifiers 95
- Malic acid
 - Chromatography on Whatman paper 12
 - Spectrophotometry 33
- Masks 110
- Master TAL 41
- Melli-mustimeter 51
- Methods for ebulliometers 22
- Micropipettes 71
- Microscopes 95
- Mini-colorimeter 31
- Mini electric compressor 17
- Mini Free Eau EVO 27
- Mini Logger USB 62
- Mustimeters 50

N-O

- Numeration cells 37
- Oxymeters 55

P

- PAL-wine 41
- Parafilm 81
- Peroxides 108
- Petri dish 37
- pH
 - buffer solutions 48
 - electrodes 17,48
 - Litmus paper 45
 - pH meter 45
 - pH tester 45
 - TITRA EVO 16
- Photometer free and total chlorine 108
- Photometer total hardness 108
- Pipette bulb 70
- Pipette controller 70
- Pipette rinsing set 70
- Pipettes
 - automatic 71
 - for musts 69
 - graduated 69
 - Pasteur 69
 - serologic 69
 - volumetric 69
- Pipette stand 70
- Pitchers 116
- Platinum electrodes 15,17
- Practical guide of alcoholometry 128

P

- Press for samples 115
- Pressure 60
- Protective faceshield 110
- Pump for pipettes 70

Q

- Quick filtration 76

R

- Rack for alcoholometers 29
- Rack for glasses 81
- RCT 0-22 41
- Reagents 121
- Refractometers
 - handheld 41
 - for bench 43
- Refrigerated baths 90
- Residual sugars
 - Fermentest 19
 - Fehling method 19
 - Spectrophotometry 33
- Ring for stand 79
- Ripper method 17
- Rods 79
- Ruler for barrel 115

S

- Safety cabinets 102
- Safety glasses 110
- Salinometers 51
- Scales 86
- SO₂
 - Extraction and determination 17,18
 - Potentiometric method 15,16
 - Ripper method 13
- Sorbic acid 12
- Spatula 81
- Spectrophotometers 34
- Spectrophotometry 33
- Spirit sampling 115
- Stainless steel temperature probes 54
- Stands 79
- Standard samples 125
- Standard solutions 120
- Standard wine 125
- Steel plates 79
- Stereomicroscope 95
- Stills
 - all degrees 23
 - plants and liquors 23
 - simple 23
- Stir bars 84
- Stoppers (grey-blue) 78
- Stoppers (screw assembly) 80
- Storage solution 48
- Sulfilyser+ 15

T

- TA01 8

Tanks

- Tartaric acid 34
- Tartaric stability 61

Test tubes

- Test tube rack 74

Thermometers

- digital 54
- enclosed scale 53
- for ebulliometer 21
- infra-red 54
- stem type 53
- thermo-hygrometer 62
- with protective sleeve 53

Thermostatic baths

- Tips for micro/macropipettes 72

TITRA EVO

- 16

Titration instruments

- Sulfilyser+ 15

- TA01, titration stand 8

- TITRA EVO 16

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- for acetic acid in vinegars 12

- for Frantz Paul 18

- for free and total SO₂, 13

- for total acidity 7

- for volatile acidity 11

Titrivin

Total acidity

- Kit for titration 7,13

- TITRA EVO 8,16

Trap for CO₂ and acid vapors

Tube water-wine for ebulliometer

Turbidimeters

- portable 56

- bench type 57

U-V

Ultrasonic baths

Vacuum filtration

Vola 2000

Volatile acidity

See also Extraction apparatus for V.A.

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Intertek Certification France, 67, boulevard
Bessières 75017 Paris - France

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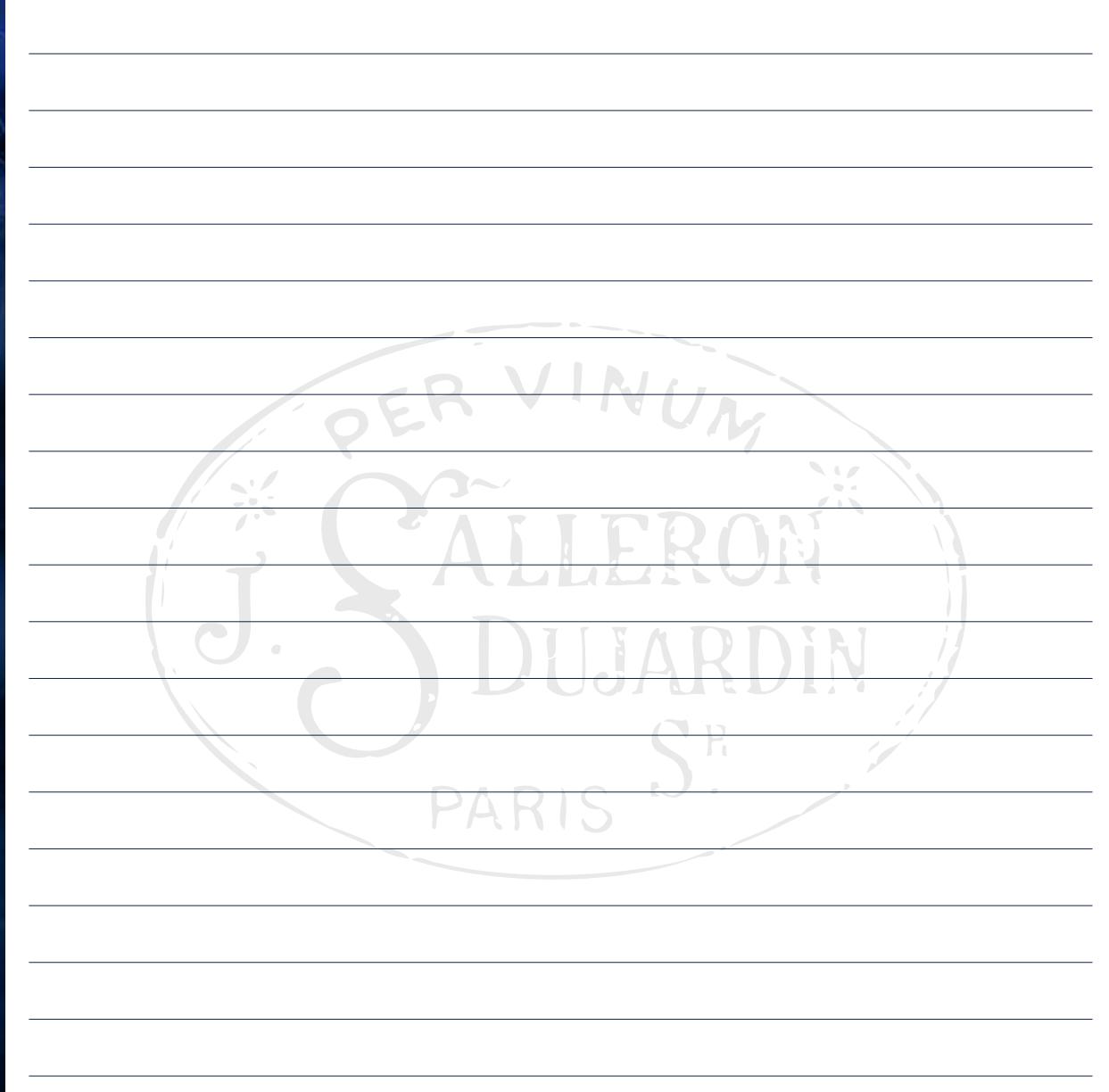
The content of this catalogue is strictly informative, and is not contractually binding for Laboratoires Dujardin-Salleron. Specifications might change at any time, without notice.
The rights of information, names and registered brands, more generally all documents on this catalogue, are reserved.
The company cannot be held responsible for errors in typography.
Photos and pictures are not contractually binding.

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© Laboratoires Dujardin-Salleron